

INSTRUCTIONS




Plastic Economy Support Stand

Catalog No. F18302-0000

To assemble the support stand,

1. Press out the small removable disc located near the edge on the top smooth surface of the base, this will create a mounting hole for the vertical rod.
2. By hand, thread one of the supplied hex nuts onto the threaded end of the stainless steel rod as far as it will go.
3. Slip the threaded end of the rod into the hole in the base and secure it by screwing the second nut onto the threaded rod protruding under the base.
4. Securely tighten the top and underside nuts with a wrench.

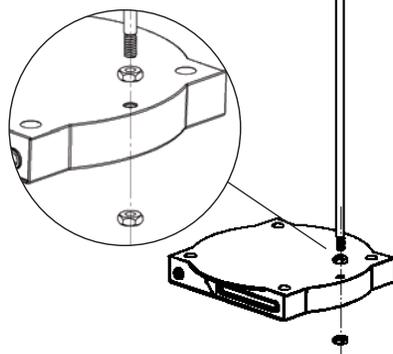

BEL-ART PRODUCTS

661 Route 23 South, Wayne, NJ 07470-6814 USA
TEL: 1-800-4BEL-ART • FAX: 973 694-7199 • www.belart.com

918302003

02/12

Scienceware® is a reg. TM of Bel-Art Products
Bel-Art Products assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products. All such advice is given and accepted at the buyer's risk.