



I N S T R U C T I O N S

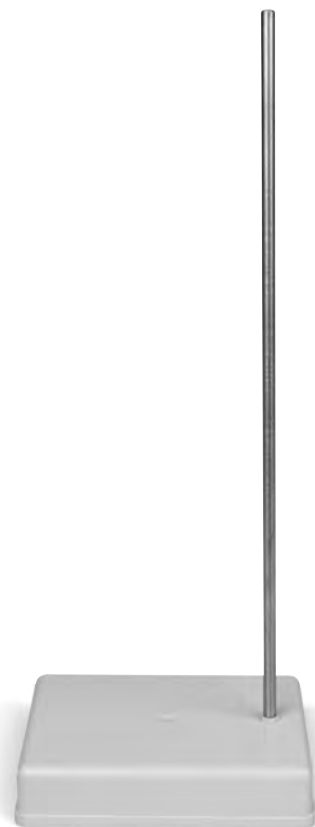

Compact Support Stand

H18303-0005

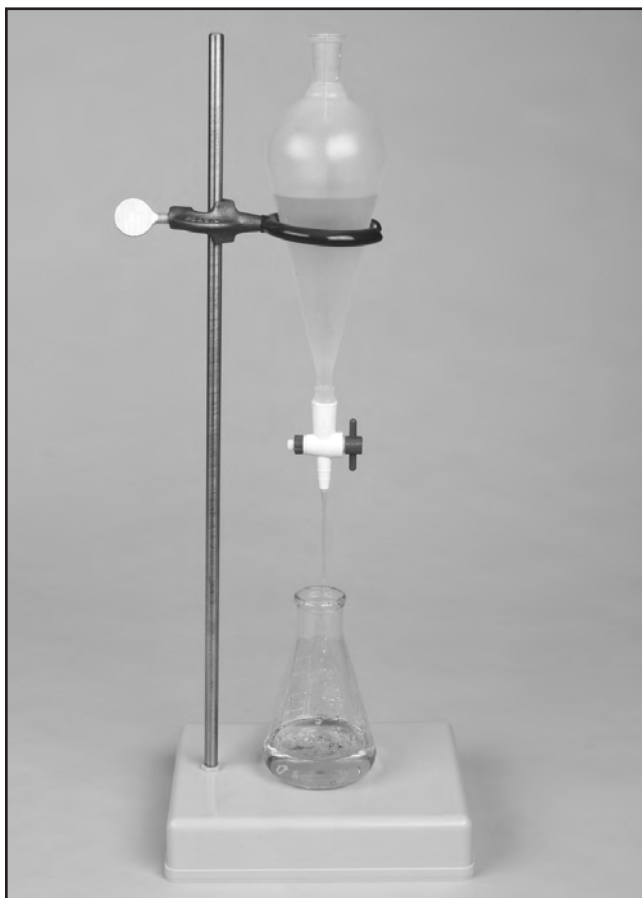

Directions for use:

1. Screw the hex nut onto the threaded end of the stainless steel rod and continue turning until it stops.
2. Screw the stainless steel rod into the desired hole in the support base and continue turning until it stops.
3. Using pliers, turn the hex nut clockwise until it makes contact with the base. Do not over tighten the hex nut as this may cause it to crack. The purpose of the hex nut is to prevent the rod from turning during use.
4. Screw the hex plug into the unused hole with a 3/16" Allen key until the top of the plug is even with the surface of the polypropylene base.

BEL-ART PRODUCTS

661 Route 23 South, Wayne, NJ 07470-6814 USA
TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • www.belart.com