

I N S T R U C T I O N S



PrepSafe™ Mini Floating MCT Rack

Inserting Microcentrifuge Tubes:

1. Hold rack securely in one hand.
2. Insert a microcentrifuge tube into the desired location.
3. Hold the tube underneath rack and slide it completely under the immobilizing tab until the tube snaps in place securing it in the rack.

Removing Microcentrifuge Tubes:


1. Hold rack securely in one hand.
2. Place your finger underneath the rack behind the tube you wish to remove and slide it out from under the immobilizing tab.

INSERTING VORTEXING ATTACHMENT:

1. Holding the rack in one hand, invert rack so that the legs are pointing upward.
2. Insert the larger end of the vortexing attachment into the hollow space within the handle of the rack until fully engaged.

REMOVING VORTEXING ATTACHMENT:

1. Holding the rack in one hand, invert rack so that the legs are pointing upward.
2. Pull vortexing attachment away from rack.

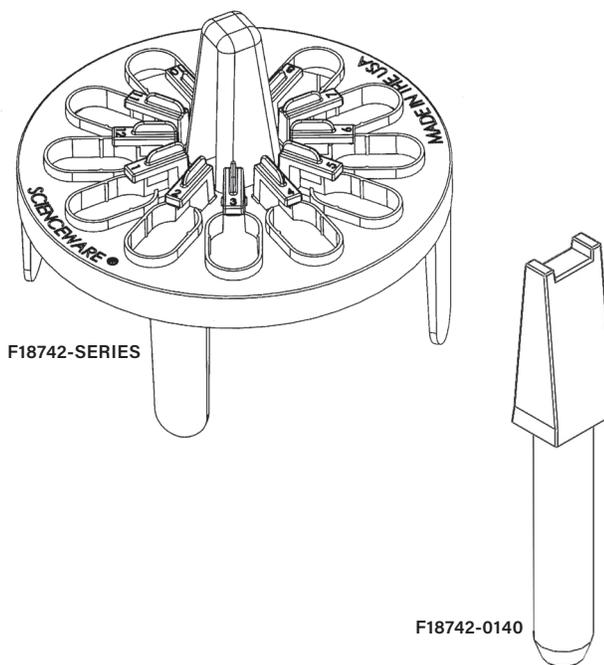

CATALOG NO.	DESCRIPTION
F18742-0120	Clear PrepSafe™ Mini MCT Rack without Vortexing Attachment
F18742-0121	Turquoise PrepSafe™ Mini MCT Rack without Vortexing Attachment
F18742-0122	Fuchsia PrepSafe™ Mini MCT Rack without Vortexing Attachment
F18742-0123	Neon Yellow PrepSafe™ Mini MCT Rack without Vortexing Attachment
F18742-0124	Clear PrepSafe™ Mini MCT Rack with Vortexing Attachment
F18742-0125	Turquoise PrepSafe™ Mini MCT Rack with Vortexing Attachment
F18742-0126	Fuchsia PrepSafe™ Mini MCT Rack with Vortexing Attachment
F18742-0127	Neon Yellow PrepSafe™ Mini MCT Rack with Vortexing Attachment
F18742-0140	Replacement Vortexing Attachment

Let others know what you think. Write a product review at www.belart.com

661 Route 23 South, Wayne, NJ 07470

TEL: 973-694-0500 • FAX: 973-694-7199 • www.belart.com

Bel-Art – SP Scienceware assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products. All such advice is given and accepted at the buyer's risk. © 2020. All Rights Reserved. Bel-Art® – SP Scienceware® is a division of SP Industries, Inc.



918742028 – Rev 2

