




No-Wire™ Half Racks

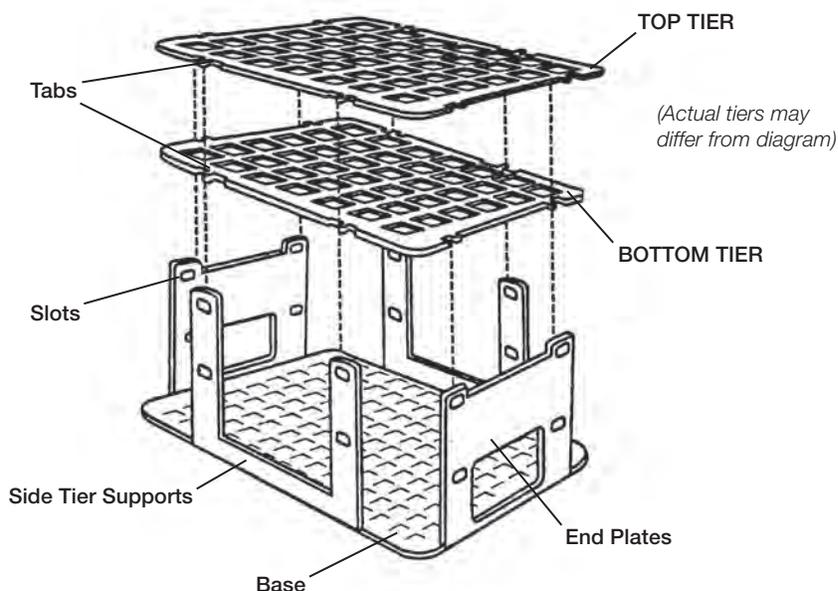


1. Place the base on a flat work surface with the word "BOTTOM" facing down.
2. Bend up one end plate. Snap the bottom tier end tabs into the lower set of slots in the end plate. Be sure the bottom tier imprinted numbers face up.
3. Position the top tier so that imprinted numbers face up and align with the numbers on the bottom tier. Snap the top tier end tabs into the upper set of slots in the end plate.
4. Bend up the opposing end plate and snap the bottom tier and then top tier end tabs into their matching slots in the end plate.

5. Bend up one side tier support and snap the bottom tier and then top tier side tabs into their matching slots in the side tier supports.
6. Bend up the remaining side tier support and snap the bottom tier and then top tier side tabs into their matching slots in the side tier supports.

Submersible and works well in water baths
Autoclavable at 250° F (121°C)
Empty racks can be stacked.
Note: Do not use with liquid nitrogen.



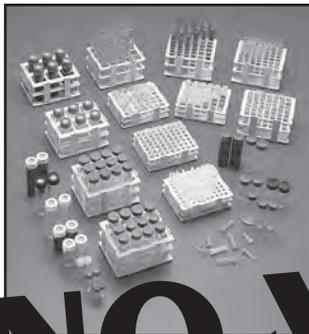

NO-WIRE™ HALF RACKS

STACKABLE • AUTOCLAVABLE • SPACE SAVER • COLOR CODED • SUBMERSIBLE
• END PLATES FOR LABELING • EASY TUBE IDENTIFICATION

NO-WIRE™ AUTOCLAVABLE VIAL AND BOTTLE RACKS

Catalog No.	Hole Size mm	Place	Rows	Fits Vials or Bottles
<input type="checkbox"/> 18514-0013	13.38	42	6 x 7	13 mm
<input type="checkbox"/> 18514-0016	16.38	30	5 x 6	16 mm
<input type="checkbox"/> 18514-0020	20.37	20	4 x 5	20 mm
<input type="checkbox"/> 18514-0025	25.40	12	3 x 4	25 mm
<input type="checkbox"/> 18514-0030	30.48	6	2 x 3	30 mm

Dimensions: 5.08" x 4.15" x 1.70" high

NO-WIRE™ AUTOCLAVABLE CUVETTE RACKS

Catalog No.	Holds Cuvettes	Place	Rows
<input type="checkbox"/> 18520-0000	10 mm	42	6 x 7

Dimensions: 5.08" x 4.15" x 1.70" high

NO-WIRE™ AUTOCLAVABLE MICRO TUBE RACKS

Catalog No.	Holds	Tubes	Place	Rows
<input type="checkbox"/> 18838-0015	1.5 ml	micro	42	6 x 7
<input type="checkbox"/> 18838-0400	.5 ml	micro	63	7 x 9
<input type="checkbox"/> 18838-0400	400µl	micro	80	8 x 10

Dimensions: 5.08" x 4.15" x 1.70" high

NO-WIRE™ AUTOCLAVABLE MICRO TUBE RACKS

Catalog No.	Holds Tubes	Place	Rows
<input type="checkbox"/> 18748-0013	13 mm	42	6 x 7
<input type="checkbox"/> 18748-0016	16 mm	30	5 x 6
<input type="checkbox"/> 18748-0020	20 mm	20	4 x 5
<input type="checkbox"/> 18748-0025	25 mm	12	3 x 4
<input type="checkbox"/> 18748-0030	30 mm	6	2 x 3

Dimensions: 5.08" x 4.15" x 2.83" high