

INSTRUCTIONS



INCUSHAKER TUBE HOLDER

Catalog No. 18798-0000

The Incu-Shaker is designed to hold various sizes of centrifuge and test tubes in a floor or bench top shaking incubator. It holds several 50 ml, 12 ml, 5 ml, 2 ml, 1.5 ml tubes [see chart]. They are positioned at an angle, which produces complete mixing/shaking for better results in any analysis. The Incu-Shaker is made of wide gauge polypropylene, suitable also for use in water baths, and is designed for a long service life in a wide variety of research applications.



Specifications:

Base Mount: 6 1/4" diameter


Tube Holder Configurations:

Tube Size	Hole Size	Number of Holes
50mL	1.40" dia.	6
12mL	0.650" dia.	12
5mL	0.465" dia.	12
2 & 1.5mL	0.420" dia.	18

Incu-Shaker Installation

The Incushaker is easily installed in your existing vibrating incubator. When you have removed the Incu-Shaker from all packing material, you will need to disassemble it in order to set it in place securely.

1. Remove each of the two plastic thumbscrews that attach the v-shaped assembly to the two support posts.
2. Place the Incu-Shaker base on the shaker mounting platform and use the four screws supplied with the Incu-Shaker to attach the base to the platform. Depending on the model of shaker, anywhere from one to four holes in the Incu-Shaker base will match up.
3. Put the v-shaped tube holder back in place. Re-install the two plastic thumbscrews.

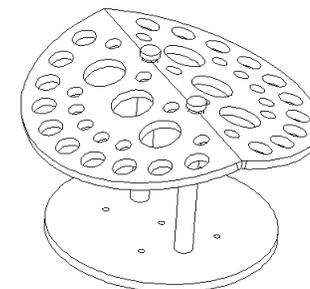

The Incu-Shaker is now fully installed.

Parts List:

The complete Incu-Shaker includes all of the following replaceable parts.

Part	Part Number	Qty. per unit
Top	91879-8101	1
Base	91879-8103	1
Post	91879-8102	2
Screw for post	99930-0302	2
Thumbscrew for post	99953-0300	2
Screw for mounting (8-32 x 1")	99905-0267	4

BEL-ART PRODUCTS

661 Route 23 South, Wayne, NJ 07470-6814 USA
TEL: 1-800-4BEL-ART • FAX: 973 694-7199 • www.belart.com

918798105

Scienceware® is a reg. TM of Bel-Art Products

Bel-Art Products assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products. All such advice is given and accepted at the buyer's risk.

03/12