

# INSTRUCTIONS




## Scienceware® Floating Rack for Cryotubes

Catalog No. F18876-0020

This 20 place round floating rack fits most cryotubes and holds them tightly in circular patterned holes, 12 in the outer ring and 8 in the inner ring. The rack floats in water and is used for quick thawing of frozen cells in 37°C water baths as well-as-for holding cryotubes during work or incubation. The filled rack is unloaded by simply pushing the rack down against the bench until the short legs touch the bench

The blue polypropylene rack is autoclavable. The rack is 1/4" (6.4 mm) thick, 3 3/4" (95mm) diameter with 3/4" (19 mm) long legs and a center handle.

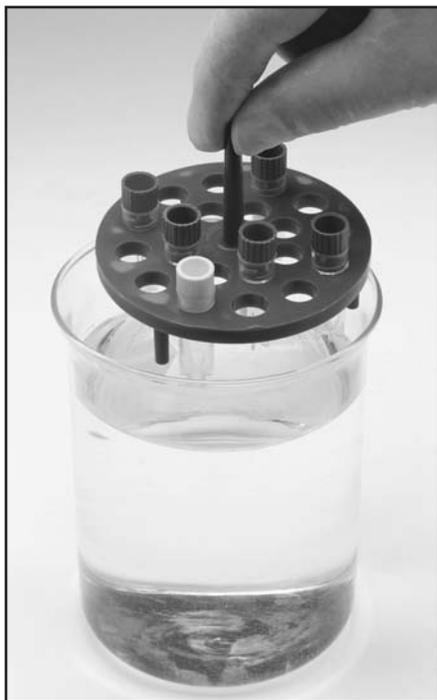

## BEL-ART PRODUCTS

661 Route 23 South, Wayne, NJ 07470-6814 USA  
TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • [www.belart.com](http://www.belart.com)

918876021

Scienceware® is a reg. TM of Bel-Art Products

Bel-Art Products assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products. All such advice is given and accepted at the buyer's risk.

03/12

*Keep Cool-Prevent Sample Loss*

## Cryo-Safe™ Cold Box

Catalog No. F18847-0002



The problem of keeping your samples cold on the benchtop or during transport between locations is solved with the new Scienceware® CryoSafe™ Cold Box.

The Scienceware® Cryo-Safe™ Cold Box is a compact unit designed to maintain a -15°C temperature for 1.5 hours (when taken out of a -20°C freezer). It will keep your samples below 0°C for four hours. A coolant is sealed into the walls of the polystyrene container. The unit is activated by placing it in a -20°C freezer for 24 hours.

The Scienceware® CryoSafe™ Cold Box has spaces for twelve 1.5mL microcentrifuge tubes. Economical and easy to use, the units are stackable to conserve freezer space.



## BEL-ART PRODUCTS

661 Route 23 South, Wayne, NJ 07470-6814 USA  
TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • [www.belart.com](http://www.belart.com)

Contact your favorite dealer or visit our web site @ [www.belart.com](http://www.belart.com) to see our complete catalog and to request your own FREE copy of our Scienceware® catalog or call 1-800-4BEL-ART.