

INSTRUCTIONS



Lab-Aire® II Wall Mounted Unit Labware Drying Rack


Catalog No. 18933-0011 • 18933-0013 • 18933-0014

WALL MOUNTING:

2 hooks or 2 #8 screws should be horizontally fastened on to a vertical mounting surface exactly 10 inches apart. In the case when screws are used, make sure the screw heads are backed off 3/16 inch from the mounting surface to allow for the thickness of the key slots of the hanging tabs.

PEG ASSEMBLIES:

On each panel, pegs should be arranged with the small pegs on the top rows and large pegs on the bottom rows. To ensure pegs are not impacted too tightly for removal or rearrangement, do not force the pegs into the peg holes.

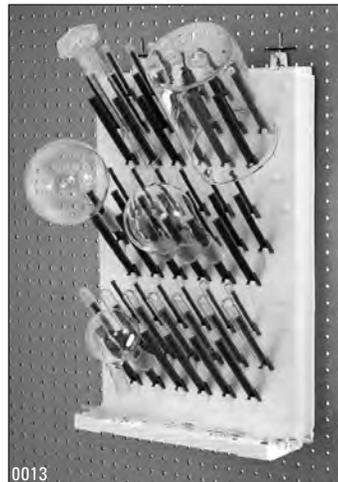

REPLACEMENT PARTS

DESCRIPTION	PART NO.
Bag of 19 Small Pegs & 6 Plugs	18931-0001
Bag of 19 Medium Pegs & 6 Plugs	18931-0002
Bag of 19 Large Pegs & 6 Plugs	18931-0003

BEL-ART PRODUCTS

661 Route 23 South, Wayne, NJ 07470-6814 USA
TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • www.belart.com

918927113

06/08

Scienceware® is a registered TM of Bel-Art Products • Bel-Art Products assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products. All such advice is given and accepted at the buyer's risk.
©2008 Bel-Art Products. All Rights Reserved.