

INSTRUCTIONS



Splash Shields

Catalog No. F24966 Series


SINGLE ARM SHIELD

1 arm with neck fitting
2 handles with bolts/washers
(these shields need 1 attachment bracket!)

TWIN ARM SHIELD

2 arms with neck fittings
2 handles with bolts/washers
(these shields need 2 attachment brackets!)

PARTS LIST/PLEXIGLAS® ASSEMBLY

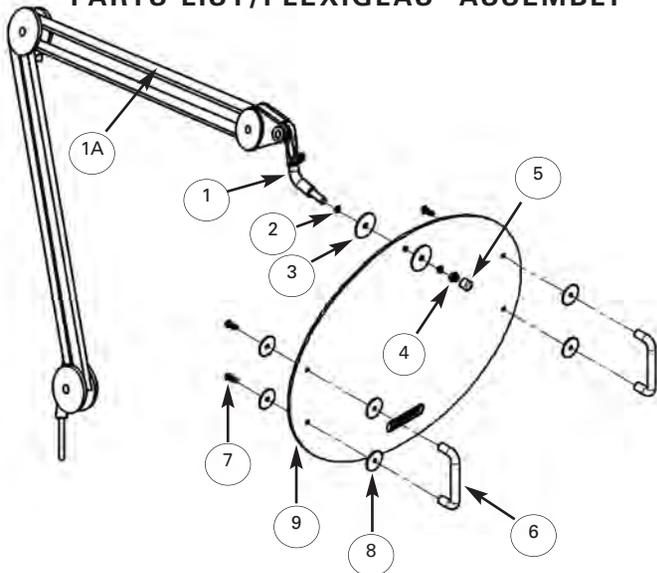

PART NUMBER AND QUANTITY

- | | |
|---------------------------|-------------------------|
| 1A. Arm (1). | 5. Cap, Red (1). |
| 1. Neck (1). | 6. Handle (2). |
| 2. Washer, Lock (2). | 7. Screw (4). |
| 3. Washer, 1.25 O.D. (2). | 8. Washer, 1" O.D. (8). |
| 4. Lock Nut (1). | 9. Shield, Oval (1). |

NOTE: ALL SHIELDS NEED ATTACHMENT BRACKETS!

WEIGHTED BASE ADAPTER
F24966-0008



See opposite side for assembly.

FIXED TABLE BRACKET
F24966-0006



C CLAMP
F24966-0005




WALL BRACKET
F24966-0007



INSTRUCTIONS

Weighted Base Bracket

Catalog No. 24966-0008

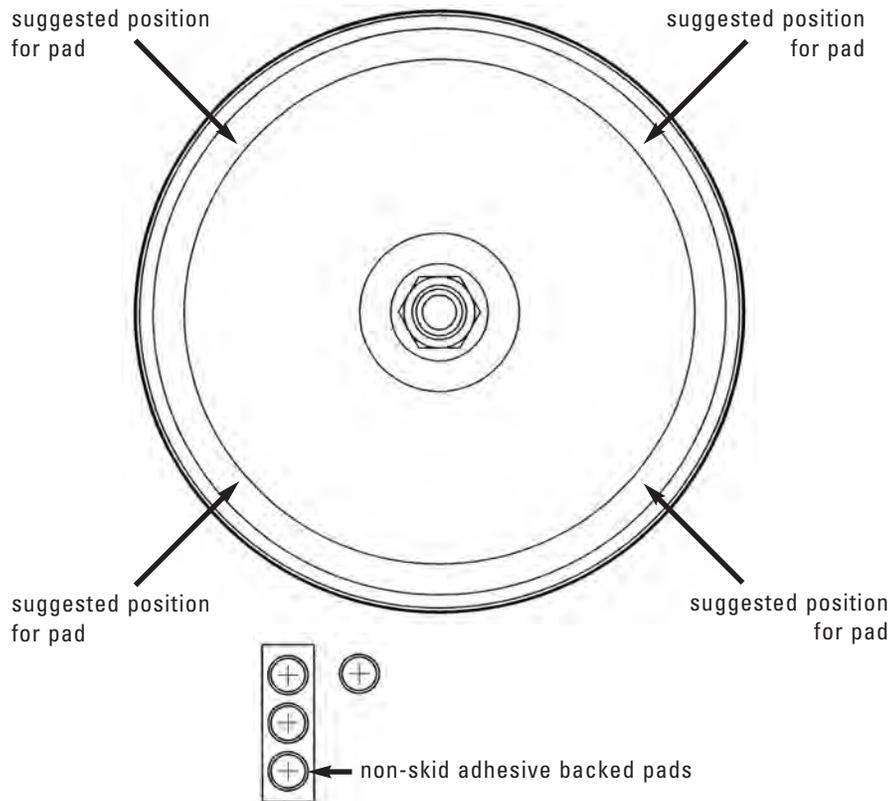

1. Apply the four (4) non-skid adhesive backed pads, included with your Scienceware® Weighted Base Bracket, to the four positions on the underside of the base.
2. Insert the shield arm into the receptor on the upper side of the weighted base.



BEL-ART PRODUCTS

661 Route 23 South, Wayne, NJ 07470-6814 USA
TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • www.belart.com

924966109

Scienceware® is a reg. TM of Bel-Art Products
Bel-Art Products assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products. All such advice is given and accepted at the buyer's risk.

08/12

