

# I N S T R U C T I O N S



## Mercury Collector Replacement Pads

Catalog No. F36251-0000

These open cell polyurethane pads are replacements for the pad supplied with the Bel-Art mercury collector F32650-0000

When handling the used pad, wear gloves and a mask to prevent contact with hazardous mercury. Dispose of the used pad in accordance with local regulations.

1. Using a scraper, such as a putty knife, remove the used pad from the three legged platform and scrape off excess dried cement.
2. To install a new pad on the platform, use rubber cement (usually available at art supply stores). Apply a layer of cement to the platform and center the pad on the platform with light pressure. For your safety, comply with the warnings on the rubber cement instructions.

661 Route 23 South, Wayne, NJ 07470



TEL: 1-800-423-5278 • FAX: 973-694-7199 • [www.belart.com](http://www.belart.com)

936251004 – 05/15

Bel-Art – SP Scienceware assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products.

All such advice is given and accepted at the buyer's risk.

© 2015. All Rights Reserved.

Bel-Art® – SP Scienceware® is a division of SP Industries, Inc.

