

INSTRUCTIONS



Bel-Bulb™ Pipettor

Catalog No. 37887-0000

The Bel-Bulb™ Pipettor fits all pipettes having top diameters of 6 to 8mm.

Hold the pipette in one hand and place the tapered chuck of the Bel-Bulb™ Pipettor over its top to create a seal. Squeeze the bulb to evacuate the air and insert the pipette tip into the liquid.

Release pressure on the bulb to draw up liquid. When the desired level is reached, quickly remove the bulb and seal the pipette with the index finger of the hand holding the pipette.

For cleaning: The chuck may be removed from the bulb and both washed in normal cleaning solutions.

NOTE: Do not autoclave.



OTHER FINE SCIENCEWARE® PRODUCTS FOR YOUR LABORATORY

PIPETTE SUPPORT RACKS

Catalog No. 18953 Series

For orderly, space-saving pipette storage. These sturdy racks holds 50 pipettes in a vertical 5 x 10 arrangement. Top and middle decks have 16mm diameter holes; bottom deck holes are tapered from 16mm to 6mm diameter. Pipette tips will not contact table surface. The end plates have handles for easy carrying. Autoclavable at 121°C (250°F). Dimensions are 213 x 114 x 222mm high (8 3/8" x 4 1/2" x 8 3/4").



F18953-0000

BEL-ART PRODUCTS

661 Route 23 South, Wayne, NJ 07470-6814 USA
TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • www.belart.com

9378876002

06/08

Scienceware®, is a registered TM of Bel-Art Products • Bel-Bulb™, is a trademark of Bel-Art Products
Bel-Art Products assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products. All such advice is given and accepted at the buyer's risk.
©2008 Bel-Art Products. All Rights Reserved.