

I N S T R U C T I O N S



Fast Release Pipette Pump™ II

Catalog Number 37911 Series

The Scienceware® Fast Release Pipette Pump™ II possesses a unique collar that requires no adjustment. The collar clamps the chuck ring to the barrel and holds the chuck in proper alignment to accept a pipette. A stabilizing mechanism inside the collar prevents the pipette from wobbling while in use.

To insert pipette:

1. Depress the plunger.
2. Insert the pipette into the collar of the pipette pump
3. Push the pipette gently past the stabilizing mechanism inside the pump's collar until it is well seated in the soft chuck. A gentle twisting motion will ease insertion of the pipette

To pipette:

1. Rotate wheel and the liquid will rise in the pipette.
2. Stop rotating the wheel when the required volume has been aspirated.
3. To release the liquid drop by drop, turn the knurled wheel slowly in the opposite direction.
4. For fast release of the liquid, press the side lever until the pipette is empty.



For Pipettes	Catalog Number
Up to 2ml	37911-1002
Up to 10ml	37911-1010
Up to 25ml	37911-1025

ATTENTION GLASS PIPETTE USERS

Caution is advised when using glass pipettes with this product. Glass may break and shatter when forced. Consult your lab safety officer for proper protocol on glass handling.

For replacement parts, please contact Bel-Art Products customer service,
(973) 694-0500

OTHER FINE SCIENCEWARE® PRODUCTS FOR YOUR LABORATORY

Write-On™ Tape Dispensers make labeling easy and fast


Producing legible labels is so much easier on a smooth, flat platform than on curved surfaces such as bottles or vials. The new Scienceware® Write-On™ Tape Dispenser provides this convenience by always having labeling tape positioned on the large writing platform, ready to make into a label. Select either the benchtop model with weighted base or economy model that can be wall mounted to save space. All models are easy to use by both left and right handed users.

Each dispenser can hold 5 or more tape rolls with widths from 12 - 25mm (½ - 1"). Ideal for use with the full line of Scienceware®, Write-On™ Tapes, F13463 series, available in 6 colors and three widths, 12, 19, and 25mm (½", ¾" and 1"). This advanced tape adheres to any clean surface and resists moisture and most solvents. It withstands long exposure to temperatures as low as -73°C (-68°F) and up to 80°C (176°F) yet peels off without leaving a sticky residue.

Contact your preferred distributor to order these and other Scienceware® products.

CATALOG NO.	DESCRIPTION	DIMENSION
F13472-0000	Benchtop Model	302 x 143 x 121 mm H (12 x 5 ⁵ / ₈ x 4 ³ / ₄ " H)
F13473-0000	Economy Model	302 x 143 x 106 mm H (12 x 5 ⁵ / ₈ x 4 ³ / ₁₆ " H)

Economy Model (shown)
Can be wall-mounted

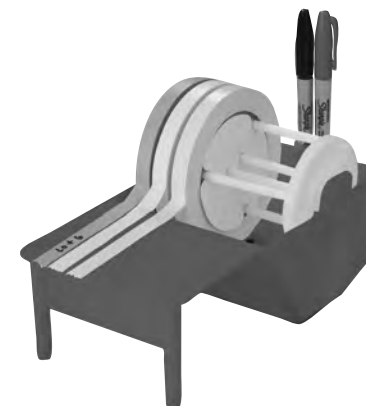

Benchtop Model (not shown)
Includes weighted base and stainless steel blade



Bel-Art Products

661 Route 23 South, Wayne, NJ 07470-6814 USA
TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • www.belart.com

937911109

02/12

Scienceware®, is a reg. TM of Bel-Art Products • Pipette Pump™ is a trademark of Bel-Art Products
Bel-Art Products assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products. All such advice is given and accepted at the buyer's risk.