



Secador® Carrying Cases Catalog No. F42070-0000 & F42070-0001

Secador Carrying Case Desiccators are molded of a durable co-polyester resin in a clear or blue color that blocks over 99% of all harmful UV light. Doors feature a patented seal and are fitted with a dial hygrometer*. A loop on the door permits a padlock or a tamper evident tag to be used to prevent unauthorized opening.

Unlike acrylic desiccators, the co-polyester resin construction of the Secador Carrying Cases can be cleaned (and disinfected) with ethanol or most laboratory detergents. Do not use abrasive cleansers as these may scratch the surface. We recommend the use of a clean, soft cloth and Bel-Art Products Aquet® Detergent, Catalog No. F17094-0030



Carrying Case should be placed on horizontal surface with the handle on top. Place desiccant cartridges under the shelf in the bottom. The desiccant will absorb moisture from the air in the case. Place items onto shelf and close the door. Take care when transporting to avoid spilling the contents.

CAUTIONS for all Secador Desiccators:

- Do not place in front of a heating or A/C duct
- Do not place near a heating device
- Do not place in areas with smoke or steam
- Do not attempt to introduce pressure or vacuum inside the cabinet
- Do not place any explosive or hazardous chemicals in the cabinets
- Do not autoclave
- Indoor use only

The user should be aware that if the equipment is used in a manner not specified by the manufacturer, the protection provided by the equipment may be impaired. All of the Secador® Desiccators are molded in sections which are then sealed tight using silicon gaskets and unexposed steel bolts. This hardware is factory assembled and no attempt should be made to alter the tightness of the nuts.

*The analog hygrometer supplied with this carry case desiccator is a reliable and accurate instrument. The hygrometer is factory calibrated. Do not attempt to adjust the settings.

BEL-ART PRODUCTS

Wayne, NJ 07470 USA • TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • www.belart.com

970001085

© 2013 Bel-Art Products. All rights reserved. Scienceware®, is a reg. TM of Bel-Art Products • Secador®, Aquet™, Cleanware™ are trademarks of Bel-Art Products

Bel-Art Products assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products. All such advice is given and accepted at the buyer's risk.

12/13

