



STERILEWARE® SENSE-ABLE SCOOPS® SAMPLING TOOLS

Thank you for your interest in Bel-Art Products Sterileware® Sense-able Scoops® Sampling Tools. FDA grade high impact polystyrene Sense-able Scoops® including fragments as small as 1.6mm³ can be detected by standard metal detection and x-ray equipment, helping you avoid quality issues and protect your product.

The enclosed sample is for testing with your x-ray/metal detection equipment. If you have any questions or need further information, please call 1-800-4BELART or email salesdept@belart.com.



Sterileware® Sense-able Scoops®

Item	Capacity	Color	Bowl Dimensions (L x W x H)	Price/Box of 100
H36902-0005	60 ml (2oz)	Black	76 x 38 x 25mm (3 x 1½ x 1")	\$157.50
H36904-0005	125 ml (4oz)	Black	91 x 46 x 31mm (3½ x 1¾ x 1¼")	241.05
H36906-0005	250 ml (8oz)	Black	127 x 51 x 51mm (5 x 2 x 2")	408.15
H36902-0006	60 ml (2oz)	Blue	7.6 x 3.8 x 2.5cm (3 x 1½ x 1")	160.65
H36904-0006	125ml (4oz)	Blue	9.1 x 4.6 x 3.1cm (3½ x 1¾ x 1¼")	246.00
H36906-0006	250ml (8oz)	Blue	12.7 x 5.1 x 5.1cm (5 x 2 x 2")	416.40


2 and 4oz scoops feature a flat bottom design and remain upright for use as weigh boats or resting on a flat surface.

METAL DETECTION REFERENCE CHART

INSTRUCTIONS ON UNDERSTANDING METAL DETECTABILITY FOR YOUR EQUIPMENT

1. On the chart below, locate the size of “sphere” used for your daily quality control sensitivity set-up.
2. Locate the area of the chart that matches your product conditions and detector settings.
3. Read the graph. Any “bar” above the “sphere” size used indicates the approximate piece size of the 4 oz. scoop that can be detected on your processing equipment.

4 oz. STERILEWARE® SENSE-ABLE SCOOPS® PERFORMANCE

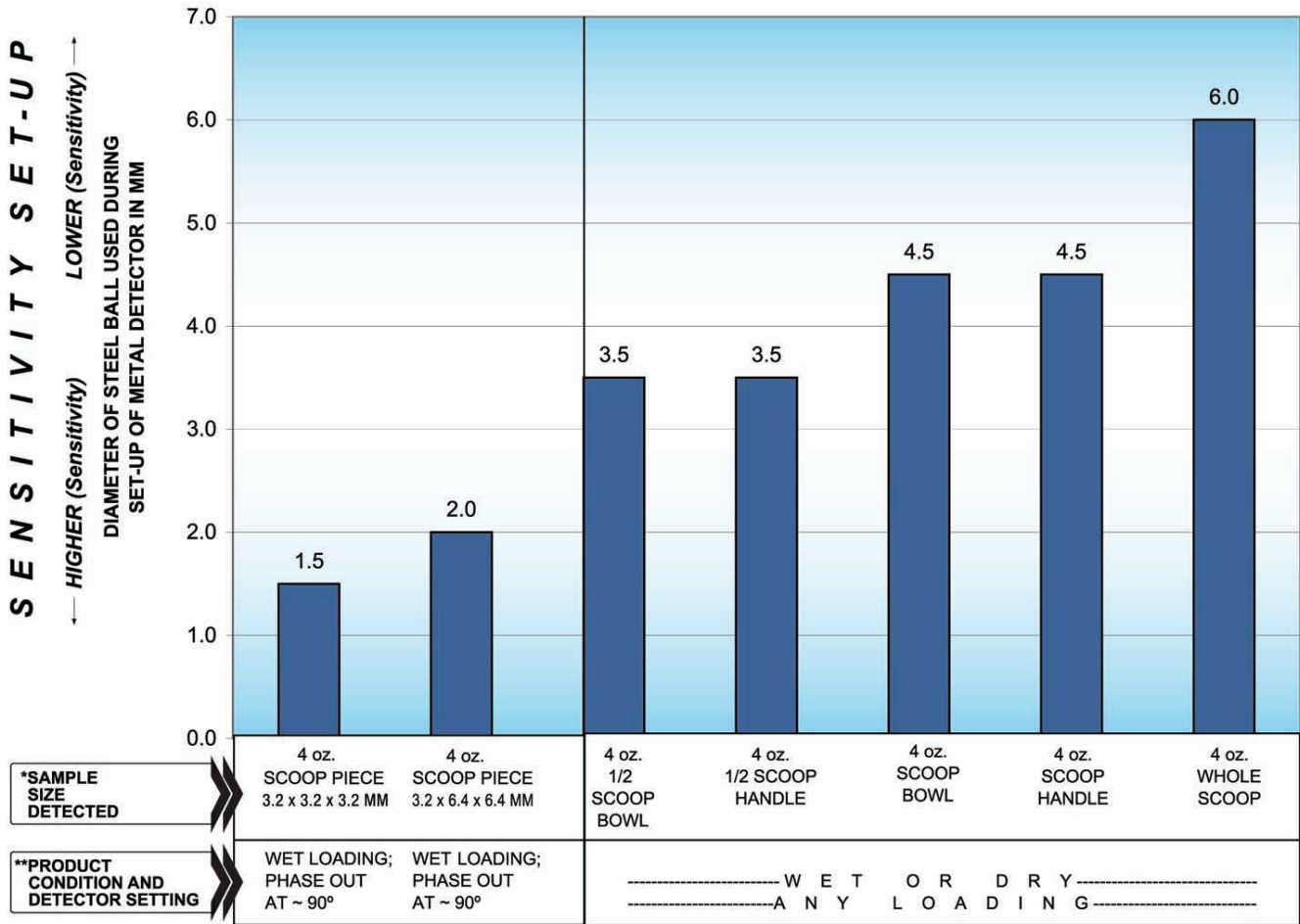

If set to higher sensitivity, the metal detector will sense a wide range of sample sizes.

All metal detection equipment and product condition(s) may differ, which may result in varying sensitivity.

* Indicates the approximate size of the sample piece that is being detected.

** Indicates your equipment setting based on the type of product that is being detected (wet or dry).

X-RAY DETECTION INFORMATION

- Scoop fragments down to 2 x 6 x 8mm are visible within an outer metal can package and even smaller fragments down to 1.6 x 1.6 x 1.6mm are visible within a plastic or glass container.
- **Results will vary and depend on the user's application.**
- **U.S. Patent No. 6,915,712 B1**

Your feedback is very important to us. In order to better serve you and improve our products, please send your feedback by email to senseablescoop@belart.com.