Instrucciones

Para ajustar la hora y calendario

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Instructions

DURAC® Multi-Lane Digital Sports Stopwatch

To set the time and calendar

Press the MODE button to display Time/Calendar
 Press the ADJUST button for 3 seconds, the seconds should blink.
 Press the START/STOP button to move from Second → Minute → Hour → Date → Month → Year to set the "blinking" digits. To adjust the digits press the LAP/SPLIT button.
 To exit the Time/Calendar set function, press the ADJUST button.
 To start measuring lap time
 Press the START STOP button
 To stop measuring lap time
 Press the START STOP button again.
 To restart measuring lap time
 Press the START STOP button

To reset stopwatch

Press the RESET button.

To start measuring multiple lap times

Press the START STOP button.
 Press the LAP/SPLIT button for first lap/split
 Press the LAP/SPLIT button for second lap/split
 Press the START STOP button to stop timing.
 Press the RESET button to reset the stopwatch.
 The time difference between the previous laps is displayed on the top most text line.

To recall Memory

Press the MODE button to view the stored lap times.
 Press either the LAP SPLIT or START STOP buttons to view the first stored lap times.
 Continue to press either the LAP SPLIT or START STOP buttons to view the next stored lap times.
 Press the ADJUST button to obtain the average of the two lap times stored.

Battery Handling/Packaging

- Remove batteries from equipment immediately after they fail to operate the equipment.
- If equipment uses two or more batteries, always replace batteries in complete sets.
- Do not attempt to recharge alkaline batteries, they are not rechargeable.
- Do not handle hot or warm batteries.
- Do not heat, incinerate, crush, puncture or mutilate batteries.
- Wear personal protective equipment if batteries show signs of leakage, bulging, swelling or deformity.

Warning: If potassium hydroxide electrolyte from an alkaline battery comes in contact with the skin, do not try to neutralize the electrolyte with vinegar or any other acidic solutions. Neutralization may trap electrolyte on the skin. Flush the affected skin area with copious amounts of water. If the battery electrolyte gets into your eyes, it can cause severe damage and/or blindness.

To register this product for a one-year limited warranty, visit www.belart.com/warranty

H-B Instrument – SP Scienceware
 Triple Accredited/Registered
 ISO 9001:2015 Registered Manufacturer
 ISO 14001:2005 Registered Manufacturer
 ISO/IEC 17025:2005 Accredited Laboratory

For more information
 973-694-0500 (p) • 973-694-7199 (f)
 info@belart.com • www.belart.com/hb

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