


INSTRUCTIONS



Vacu-Rack

Catalog No. 18977-0000

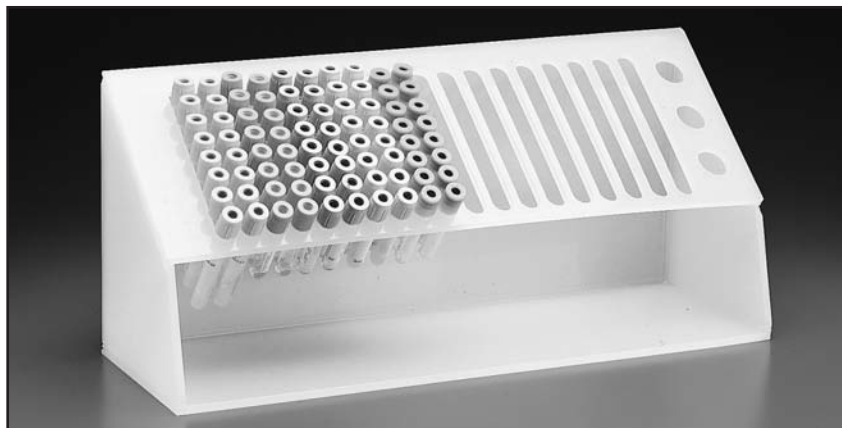

Made of polypropylene the Vacu-Rack is autoclavable and can be cleaned by chemical means.

The Vacu-Rack keeps vacuum type syringe tubes in neat order and readily available. Hang the tubes in the rack filling each of the channels. Remove the lowest tube and those behind it will slide down to keep the first row available. The tubes hang on their colored stoppers which are always visible.

The rack has slots for 1/2" (12.7mm) diameter tubes and 5/8" (15.9mm) diameter tubes. The three side holes are for needle holders.

BEL-ART PRODUCTS

661 Route 23 South, Wayne, NJ 07470-6814 USA
TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • www.belart.com

999320831

03/12

Scienceware®, is a reg. TM of Bel-Art Products • Cryo-Safe™ is a trademark of Bel-Art Products
Bel-Art Products assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products.
All such advice is given and accepted at the buyer's risk.

OTHER FINE SCIENCEWARE® PRODUCTS FOR YOUR LABORATORY

Cryo-Safe™ Cold Box

Catalog No. 18847-0002



Keep Your Microtubes Cool

The Cryo-Safe™ Cold Box provides a convenient means of preventing expensive enzymes and critical samples from warming. This handy unit consists of a molded polystyrene container with cover. There is an internal 12 place support for 1.5 ml microcentrifuge tubes. The space surrounding the tubes is filled with a sealed-in insulation material that maintains the low temperature required for safe storage and transport. When the container is taken out of a -20°C freezer it will maintain a -15°C temperature for 1.5 hours and will maintain a temperature below 0° C for four hours.



BEL-ART PRODUCTS

661 Route 23 South, Wayne, NJ 07470-6814 USA
TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • www.belart.com