

INSTRUCTIONS




Adapter and Trigger Sprayer Catalog No. F11620-0050

The Bel-Art – SP Scienceware® Adapter has one end to fit the opening on all 500ml and 1000ml Bel-Art – SP Scienceware® wash bottles and another end to fit the Bel-Art – SP Scienceware® Trigger Sprayer. The adapter and trigger sprayer are compatible with a wide range of chemicals. However, because the sprayer can be adjusted to produce a fine mist, it is not recommended for use near open flames or other sources of combustion. Directions for affixing the adapter and trigger sprayer are provided below.

Directions for Assembly:

1. To attach adapter to a 1000ml Bel-Art – SP Scienceware® wash bottle, skip to Step #7
2. To attach adapter to a 500ml Bel-Art – SP Scienceware® wash bottle, you will need to cut the tubing as described below
3. Grasp the tubing and pull it away from the trigger sprayer to separate it from trigger sprayer
4. Align filter end of the tubing with line drawing shown in Figure 2
5. Using scissors, cut tubing to the length corresponding with Figure 2
6. Re-insert tubing into trigger sprayer
7. Screw adapter onto bottle
8. Guide the tubing of the trigger sprayer through the opening on the adapter and into the bottle
9. Screw trigger sprayer into adapter
10. The spray adjustment pattern can be varied from mist to stream by turning the spray adjustment nozzle as shown in Figure 1

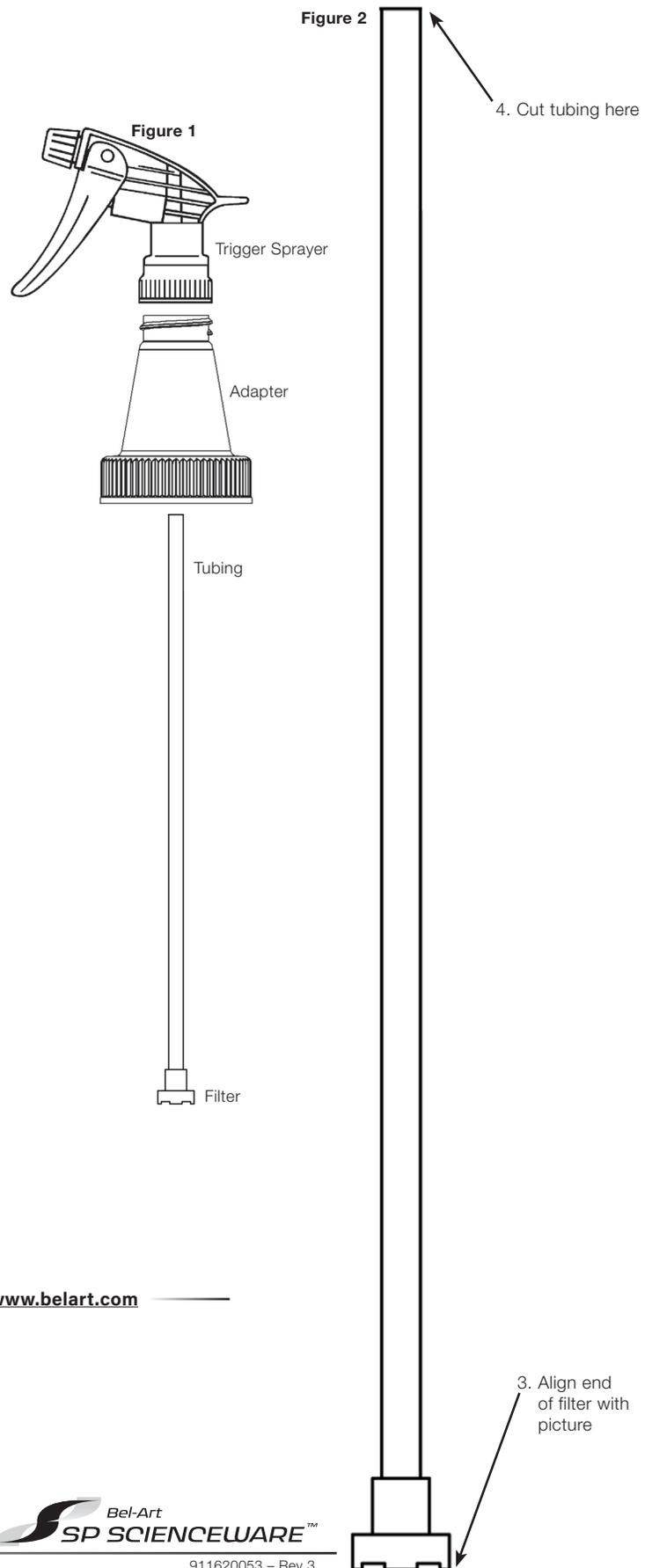

Let others know what you think. Write a product review at www.belart.com

661 Route 23 South, Wayne, NJ 07470

TEL: 973-694-0500 • FAX: 973-694-7199 • www.belart.com

Bel-Art – SP Scienceware assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products.

All such advice is given and accepted at the buyer's risk.

© 2019. All Rights Reserved.

Bel-Art® – SP Scienceware® is a division of SP Industries, Inc.



911620053 – Rev 3

