

INSTRUCTIONS





Clavies® Biohazard Bag Holders

Catalog No. F13192-0001, -0002, -0003

1. Small F13192-0001 (7.3 x 14 x 18.5")

Parts List

1. Tray (1)
2. Wire Frames (2)
3. Plastic Wire Anchors (4)

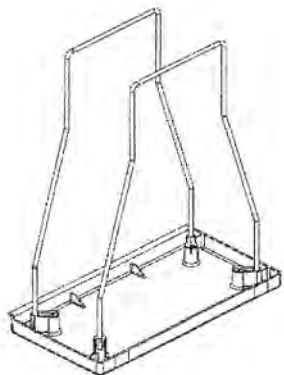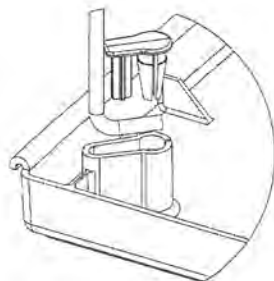

- Turn the tray (Part #1) right side up so the four wells are visible.
- Push the ends of the stainless steel wire frames (Part #2) into the wells.
- Push the four plastic wire anchors (Part #3) into the wells until they are fully seated on the rim of the well and lock the wire frame into place.
- Drape the appropriate size bag over the wire frames and roll the edge over the top of the frames.

Recommended bags for small Clavies® Biohazard Bag Holder are Scienceware® F13160-0009, F13160-0005, H13185-1224 & F13164-1923

BEL-ART PRODUCTS

661 Route 23 South, Wayne, NJ 07470-6814 USA
TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • www.belart.com

913192011


07/09

Scienceware® and Clavies® are reg. TM of Bel-Art Products
Bel-Art Products assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products. All such advice is given and accepted at the buyer's risk.


2. Medium F13192-0002 (14 x 14 x 21.5"), Large 13192-0003 (14 x 14 x 27.2")

Parts List

1. Trays (2)
2. Wire Frames (2)
3. Plastic Wire Anchors (8)
4. Wire Locking Clips (2)

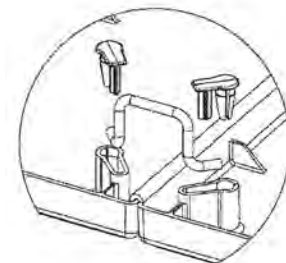

- Turn the two trays (Part #1) right side up so the eight wells are visible.
- Join the two trays together by slipping the rolled edge of one over the straight edge of the other.
- Install the two stainless steel wire locking clips (Part #4) in the center well to lock the trays together.
- Push the ends of the stainless steel wire frames (Part #2) into the wells.
- Push the eight plastic wire anchors (Part #3) into the eight wells until they are fully seated on the rim of the well and lock the wire frames and wire locking clips into place.
- Drape the appropriate size bag over the wire frames and roll the edge over the top of the frames.

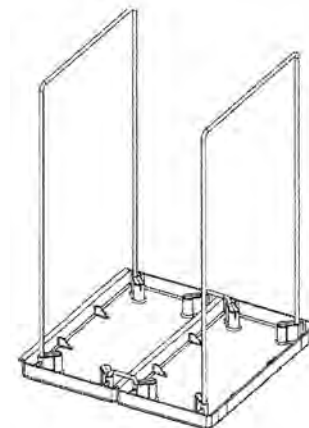

*Recommended bags for medium Clavies® Biohazard Bag Holder are Scienceware® F13161-0009, F13161-0005 & H13185-2430
Larger bags can also be substituted.*

*Recommended bags for Large Clavies® Biohazard Bag Holder are Scienceware® F13162-0009, F13162-0005, H13185-2436, F13164-2535, F13165-2535, F13164-3138 & F13165-3138
Larger bags can also be substituted.*



**Improve sanitation by adding a cover to the bag holder.
Recommended cover for the medium or large Clavies®
Biohazard Bag Holder is F13192-0102.**

**For more information on this item, please visit www.belart.com
and search for this item number**

