



Touch-Free™ Automatic Waste Can

Catalog No. F13202-0010 (2.5 gallon), F13202-0020, F13202-0022 (7.3 gallon) & F13202-0030 (12 gallon)

The patented Touch-Free™ Automatic Waste Can opens automatically when your hand or object approaches within 8 inches of the built-in sensor in the lid and closes automatically after 3 seconds. It is safe and simple to operate - ideal for laboratories, hospitals, physician and other medical offices.

Included with each unit are 4 labels: Biohazard, Recycle, Trash and a blank "Label Your Own" label. Select the label that is appropriate for your use and place it on the lid to clearly identify the purpose of the waste can and prevent safety and disposal problems caused by mixed waste.

Scienceware® Biohazard Disposal Bags, size 14" x 19", fit perfectly in the 2.5 gallon model. Ask for part number F13164-1419 or F13165-1419 (for extra strength model).

Scienceware® Biohazard Disposal Bags, Clavies® Autoclavable Bags and Bag Liners measuring 24 x 30", 24 x 36" or 25 x 35" fit will fit 7.3 and 12 gallon models.

Size Catalog No.

- 24 x 30" F13161-0009 or F13161-0005
(Biohazard bags with or without sterilization indicator patch)
- 24 x 36" F13162-0009 or F13162-0005
(Biohazard bags with or without sterilization indicator patch)
- 25 x 35" F13164-2535 or F13165-2535
(Standard or extra strength Biohazard bags)
- 24 x 30" H13185-2430 (Clavies® Autoclavable Bags)
- 24 x 36" H13185-2436 (Clavies® Autoclavable Bags)

Set Up

- 1) Before assembly, locate power switch on the rear side of the top and turn to OFF (marked "O").
- 2) Battery compartment is located under the lid. Insert "D" size batteries (not included) and close the battery case.
- 3) Place the bag liner inside pail.
- 4) For the 7.3 & 12 gallon models, place the top onto the pail. The

top should be positioned on the pail in such a way that the power switch is located over the side of the pail that is slightly wider. Your Touch-Free™ Automatic Waste Can is now assembled and ready for use.

Operation

- 5) Press the power switch to ON (marked as "I"). The red indicator light on the lid will illuminate after 4 seconds. It will then flash every 6 seconds to indicate proper functioning.
- 6) To use the Touch-Free™ Automatic Waste Can, move your hand toward the front of the lid. The light will turn green and the lid will open automatically. The lid will remain open until your hand moves away. The lid then closes automatically after 3 seconds.
- 7) Keep Open Feature: Press the "Open" button on the left front part of the lid to keep the lid open for extended time periods. Press the "Close" button on the right front part of the lid to disable this feature and return to automatic mode.
- 8) Battery Indicator: When the red flashing light changes to an amber color, the batteries should be changed. The battery compartment is located under the lid.

Operation

- 9) Turn off the power if the Touch-Free™ Automatic Waste Can will be idle for long periods of time.
- 10) Do not open or close the lid manually. Instead, use the "OPEN" or "CLOSE" buttons, if necessary.
- 11) Use cleaning cloths or wipes to clean the removable top. Turn unit OFF and remove batteries BEFORE any cleaning operations are performed. The pail can be washed and dried separately as needed.

661 Route 23 South, Wayne, NJ 07470

TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • www.belart.com

Scienceware® is a registered TM of Bel-Art Products.
Bel-Art Products assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products.
All such advice is given and accepted at the buyer's risk.
© 2013 Bel-Art Products. All Rights Reserved.



913202019 - 07/13

