



## Weighted Beaker/Flask Holders Catalog No. F18309-0100, 0250, 0600, 1000

Scienceware® Weighted Beaker/Flask Holders grip beakers or flasks for immersion in water baths. Circulating water in water baths may cause the beaker or flask to tip or drift; partially filled vessels may float. Use of the weighted holders prevents such accidental dislodging.

1. The straight prongs on the bottom of the stainless steel clamp are inserted into the center hole of the vinyl-coated lead ring.
2. Bend the prongs around the weight so the weight cannot fall off.
3. Select the size of the holder to fit the beaker or flask you are using.
4. Press down gently on the glass or plastic beaker, or flask, to insert it into the stainless steel clamp. Take normal precautions in handling glass to avoid breakage and injury.
5. Place the filled vessel with weighted holder into the water bath.
6. To release the vessel from the holder, hold down on the weight and pull up on the vessel.



## OTHER FINE SCIENCEWARE® PRODUCTS FOR YOUR LABORATORY

### Vikem® Vinyl Coated Lead Wire

*Stabilizes Vessels In A Water Bath*

This lead wire is flexible and conforms to any shape. It can be coiled into a helix, spiral, cone or countless other configurations. Curl around flasks, thread through handles or loop around bottle necks. The lead wire is 1/4" (6.4mm) in diameter and is coated with a chemical-resistant red Vikem® vinyl. The wire's flexibility makes it a very versatile weight. Length 12" (300mm), weight approximately 1/4lb. (114gm).

**18324-0000**



For more information, visit [www.belart.com](http://www.belart.com)

Help others choose the right product. Write a review at [www.belart.com](http://www.belart.com)

661 Route 23 South, Wayne, NJ 07470

TEL: 973-694-0500 • FAX: 973-694-7199 • [www.belart.com](http://www.belart.com)

Bel-Art – SP Scienceware assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products. All such advice is given and accepted at the buyer's risk. © 2020. All Rights Reserved.

Bel-Art® – SP Scienceware® is a division of SP Industries, Inc.



918309004 – Rev 2

