

# I N S T R U C T I O N S



## 15 and 50ml Tube Holder

F18796-0000



### Initial setup:

1. Use a twisting action or scissors to remove the tube supports from the base at the points outlined in the above picture,
2. Insert tube supports into the slots on the base to produce either a 15 or a 50ml tube holder.

### Directions for use:

1. Place tube between supports, ensuring that conical portion of tube is sitting in the hole in the center of the base.

**Note:** The two inner slots accommodate a 15ml tube and the two outer slots accommodate a 50ml tube.

2. Apply slight pressure to tube supports as shown in the top right photo.
3. Keep slight pressure applied to the tube supports while transporting the tube within the holder.

## BEL-ART PRODUCTS

661 Route 23 South, Wayne, NJ 07470-6814 USA  
TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • [www.belart.com](http://www.belart.com)