

# I N S T R U C T I O N S



## ADJUSTABLE ELUTION RACK

Catalog No.	Places	Tube Size
F18882-0001	42	13mm
F18882-0002	30	16mm
F18882-0003	12	25mm
F18882-0004	6	30mm



### RACK SET UP:

1. Lift front of half rack and pull out of the unit.
2. Insert the collection tubes.
3. Reinsert half rack into unit.
4. Insert elution columns into grid.
5. Adjust the height of the rack by turning the height adjustment wheel located on the handle.
6. Once the desired rack height has been reached, secure the rack height by hand-tightening the two plastic screws located on the left and right side of the rack handle.
7. Steam Autoclavable at 121°C (250°F).

### CONVERSION KITS



*Includes half rack and two grids*

Catalog No.	Places	Size
<b>A.</b> F18882-0005	42	13mm
<b>B.</b> F18882-0006	30	16mm
<b>C.</b> F18882-0007	12	25mm
<b>D.</b> F18882-0008	6	30mm

### REPLACEMENT GRIDS

*Set of two grids*

Catalog No.	Places	Size
<b>A.</b> 91888-2013	42	13mm
<b>B.</b> 91888-2016	30	16mm
<b>C.</b> 91888-2025	12	25mm
<b>D.</b> 91888-2030	6	30mm

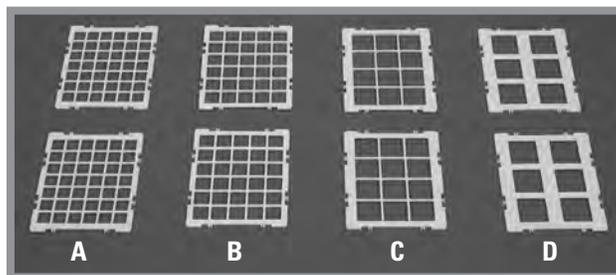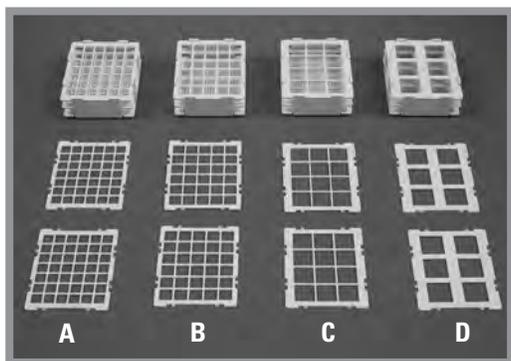

## BEL-ART PRODUCTS


661 Route 23 South, Wayne, NJ 07470-6814 USA  
 TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • www.belart.com

# I N S T R U C T I O N S





## ADJUSTABLE ELUTION RACK CONVERSION KIT

Catalog No.	Places	Size
<b>A.</b> F18882-0005	42	13mm
<b>B.</b> F18882-0006	30	16mm
<b>C.</b> F18882-0007	12	25mm
<b>D.</b> F18882-0008	6	30mm

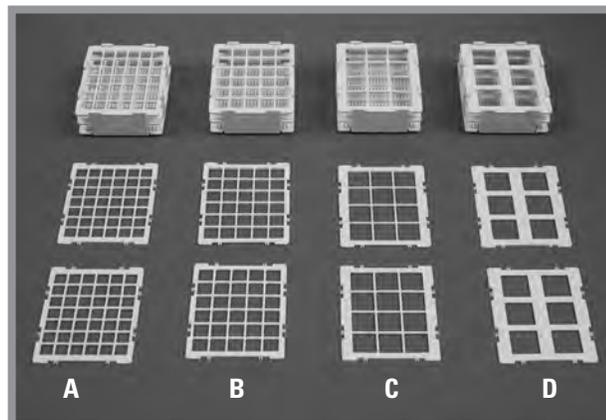

1. Lift front of the half rack and pull out of the unit. Insert new half rack at a slight angle.
2. Insert a grid into the track located immediately above the half rack.
3. Insert the second grid on the track at the very top of the rack by completing the following steps:
  - 3a. Hold the top grid in place while unfastening the two hinges.
  - 3b. Remove the old grid. Insert the replacement grid into the track. Hold the new grid in place and close the two hinges until they snap onto locking pins.

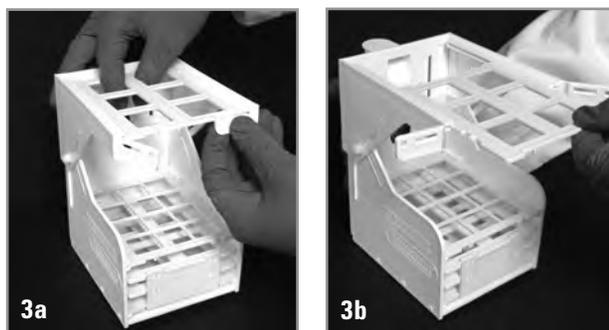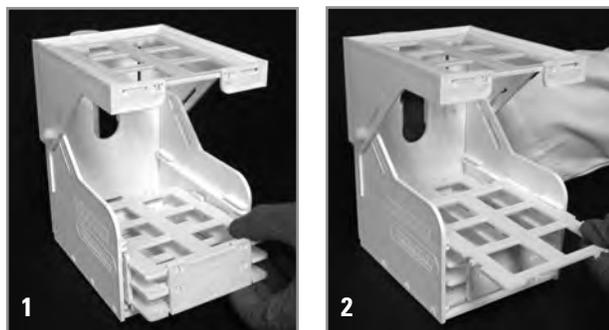

### BEL-ART PRODUCTS


661 Route 23 South, Wayne, NJ 07470-6814 USA  
TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • [www.belart.com](http://www.belart.com)

# I N S T R U C T I O N S



## ADJUSTABLE ELUTION RACK REPLACEMENT GRIDS

Part No.	Places	Tube Size
<b>A.</b> 91888-2013	42	13mm
<b>B.</b> 91888-2016	30	16mm
<b>C.</b> 91888-2025	12	25mm
<b>D.</b> 91888-2030	6	30mm

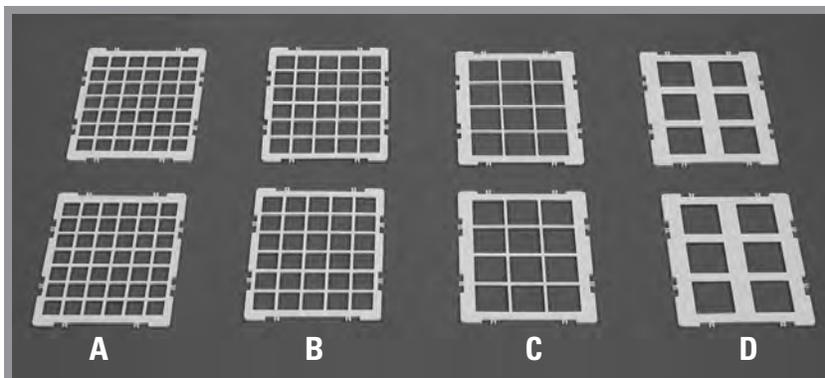

### RACK SET UP:

1. Insert a grid into the track located immediately above the Half Rack.
2. Insert the second grid on the track located on the very top of the rack by completing the following steps:
  - 2a. Hold the top grid in place while unfastening the two hinges.
  - 2b. Remove the old grid. Insert the replacement grid into the track. Hold the new grid in place and close the two hinges until they snap into place.



## BEL-ART PRODUCTS

661 Route 23 South, Wayne, NJ 07470-6814 USA  
TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • [www.belart.com](http://www.belart.com)