

I N S T R U C T I O N S

Bel-Art® – SP Scienceware® Magnetic Bead Separation Racks

WARNING!

These separation racks contain powerful Neodymium magnets. Misuse can result in injury.

- Do not use this product if you have a pacemaker or any other medical magnetized implant
- Keep away from tools and objects that could be damaged by the magnetic field
- Racks should be handled and stored away from devices subject to magnetic damage including credit cards and other products containing magnetic recording devices
- Keep away from delicate instruments, watches, electronic equipment, displays and monitors
- Magnets may attract steel or other magnetic material with high mechanical forces
- Avoid contact between two magnets
- Do not pull the magnets apart if contact has been made; twist off to prevent damage to the unit or fingers

Cleaning and Disinfection:

Should the surfaces of your SP Scienceware acrylic products need cleaning, **DO NOT USE ANY SOLVENTS** (alcohol, acetone, etc..) or abrasives. Use of these materials **will cause cracking and/or crazing of the acrylic**. For general cleaning, wash the surfaces with a mild detergent and soft cloth, then rinse thoroughly with water.

IMPORTANT — Do not autoclave

Magnetic Bead Separation Rack
for 50ml Conical Tubes
catalog number #F19900-0000




Magnetic Bead Separation Rack
for 5-15ml Conical Tubes
catalog number #F19900-0001




Magnetic Bead Separation Rack
for 1.5-2ml Centrifuge Tubes
catalog number #F19900-0002



Magnetic Bead Separation Rack
for 96-well Conical-Bottomed Plates
catalog number #F19900-0003

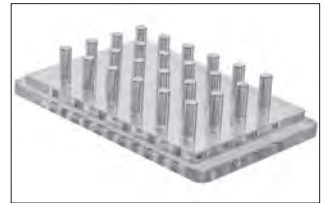

Magnetic Bead Separation Rack
for 96-well Flat-Bottomed Plates
catalog number #F19900-0004

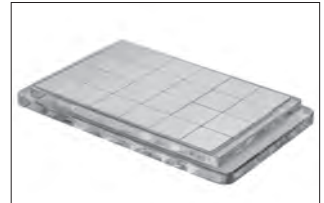

You May Also be Interested in: Bel-Art® – SP Scienceware® Micro Tube Homogenizer System

Micro Tube Homogenizer – Includes Cordless Motor,
2 AA Batteries and Pestle Adapter


F65000-0000 Cordless Homogenizer Unit, –
includes 2 AA batteries

Individual Pack

Catalog No.	Description	Case Qty.
F65000-0001	0.5 mL Pestle	100
F65000-0003	0.5 mL Tube	100
F65000-0005	0.5 mL Pestle and Tube	100
F65000-0002	1.5 mL Pestle	100
F65000-0004	1.5 mL Tube	100
F65000-0006	1.5 mL Pestle and Tube	100

Bulk Pack

Catalog No.	Description	Case Qty.
F65000-9001	0.5 mL Pestle	100
F65000-9003	0.5 mL Tube	100
F65000-9005	0.5 mL Pestle and Tube	100
F65000-9002	1.5 mL Pestle	100
F65000-9004	1.5 mL Tube	100
F65000-9006	1.5 mL Pestle and Tube	100

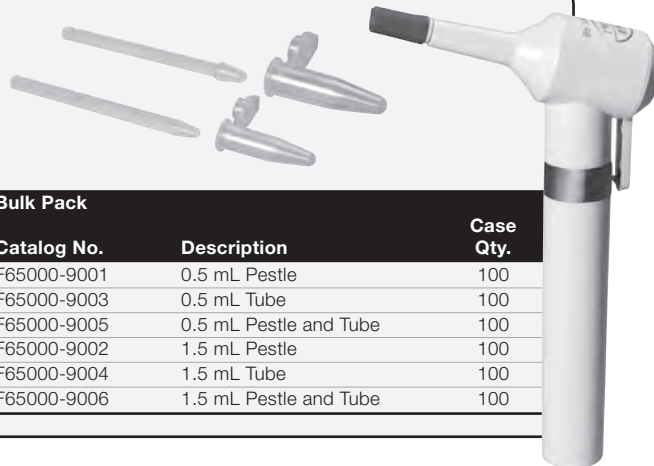

661 Route 23 South, Wayne, NJ 07470

TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • www.belart.com

Bel-Art – SP Scienceware assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products.

All such advice is given and accepted at the buyer's risk.

© 2015. All Rights Reserved.

Bel-Art® – SP Scienceware® is a division of SP Industries, Inc.

Bel-Art • H-B Instrument
SP SCIENCEWARE™

919900023 – 02/15

