

# I N S T R U C T I O N S



## Maintenance of Bel-Art – Scienceware® Eye Wash Bottles

Catalog No. F24850-0000, F24851-0000

1. Wash the Eye Wash Bottle and all accessory parts: cap, tubes, etc. in mild detergent solution.
2. Rinse well to remove all traces of detergent.
3. Under clean conditions, fill Eye Wash Bottle completely.
4. Replace top, close cap and return to emergency station.
5. If an eye wash bottle is used, empty it completely and proceed through steps 1-4 before replacing in station.
6. After using an eye wash bottle, the injured individual should see a physician at once for examination and treatment.

To ensure cleanliness, it is advisable to establish a regular eye wash bottle cleaning and eye wash solution replacement procedure. See the refilling guidelines below.



## Guidelines for Refilling Bel-Art – Scienceware® Eye Wash Bottles

The below recommendations for solution replacement are based on information gathered from sources outside Bel-Art – SP Scienceware to help assist our customers. Bel-Art Products does not guarantee the accuracy of the information gathered.

Rinse solution and frequency of change:

- *Water only - change once a week.*
- *Water and Preservative (varies by manufacturer) - change every two months.*
- *Pre-mixed Eye Wash Solution - change every 4 to 6 months.*

## Eye Wash Station and Refill



**Catalog No. F24868-0000**

### Bright, Easy to Locate

Our easily recognized Eye Wash Stations come complete with eye wash bottles, mirror, and instructions for use and service. A bracket holds the bottles securely in place, yet the bottles are easily released when pulled out. Includes station and empty eye wash bottles. We recommend filling with sterile eyewash solution catalog number 24879-9032.

**Catalog No.**

- F24865-0000
- F24866-0000
- F24868-0000

**Bottle Size**

- Station with one 500ml (16oz) empty eye wash bottle
- Station with one 1000ml (32oz) empty eye wash bottle
- Station with two 1000ml (32oz) empty eye wash bottles

661 Route 23 South, Wayne, NJ 07470

TEL: 973-694-0500 • FAX: 973-694-7199 • [www.belart.com](http://www.belart.com)

Bel-Art – SP Scienceware assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products. All such advice is given and accepted at the buyer's risk.

© 2019. All Rights Reserved.

Bel-Art® – SP Scienceware® is a division of SP Industries, Inc.



924851013 – Rev 3

