


# INSTRUCTIONS



## Liquid Sampler

Catalog Number 37827-0000

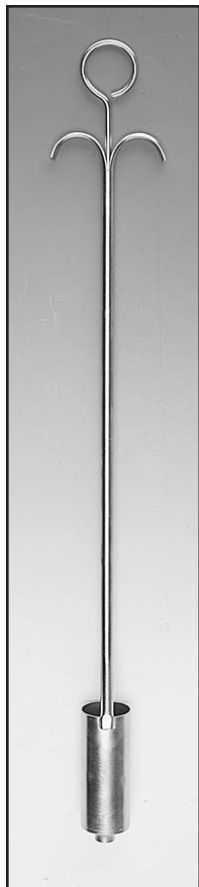

1. Be sure that the sampler is clean and free of foreign matter that might be left from packing materials.
2. The Sampler has 2 finger grips and a center ring.
3. Hold the finger grips with two fingers and put the thumb through the center ring.
4. Squeeze the center ring downward to make sure that the plunger closes tightly at the bottom.
5. Insert the cup of the Sampler into the liquid.
6. Withdraw the Sampler from the fluid and hold it over a beaker or other collecting vessel.
7. Raise the ring with the thumb to release the fluid.
8. The plunger may be unscrewed from the center rod for easier cleaning.

## BEL-ART PRODUCTS

661 Route 23 South, Wayne, NJ 07470-6814 USA  
TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • www.belart.com

937827911

03/02

Scienceware®, is a reg. TM of Bel-Art Products  
Bel-Art Products assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products. All such advice is given and accepted at the buyer's risk.

## OTHER FINE SCIENCEWARE® PRODUCTS FOR YOUR LABORATORY

### Long-Handled Dippers



For easy sampling from tanks, vats, lakes, etc., these high-density polyethylene dippers will not rust or corrode and provide excellent chemical resistance. 29mm (1 1/8") diameter handle is welded to the rigid bowl at a 45 degree angle. The 3.7M (12 ft) handle is available in one piece or two-piece screw-together construction. The end of the handle has a hook for hanging. Orders may be customized by substituting scoops in place of cups. Other styles made to order. 12 per case.

<u>CATALOG NO.</u>	<u>BOWL SIZE</u>	<u>HANDLE LENGTH</u>	<u>HANDLE</u>
F36780-0016	500cc (16oz)	91cm (3 ft)	1 piece
F36781-0016	500cc (16oz)	183cm (6 ft)	1 piece
F36782-0016	500cc (16oz)	366cm (12 ft)	2 pieces
F36782-0116	500cc (16oz)	366cm (12 ft)	1 piece*
F36780-0032	1000cc (32oz)	91cm (3 ft)	1 piece
F36781-0032	1000cc (32oz)	183cm (6 ft)	1 piece
F36782-0032	1000cc (32oz)	366cm (12 ft)	2 pieces
F36782-0132	1000cc (32oz)	366cm (12 ft)	1 piece*

\*One piece handles measuring 12 ft. cannot be shipped via small package carrier.



## BEL-ART PRODUCTS

661 Route 23 South, Wayne, NJ 07470-6814 USA  
TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • www.belart.com