

INSTRUCTIONS



Mini-Vacuum Desiccator

Catalog No. 41099-0000

The Mini-Vacuum Desiccator consists of a clear top section, a white bottom section, an O-ring and a vacuum retention valve.

The vacuum retention valve fits on the tapered neck at the top. It accepts 1/4" inside diameter tubing.

Place the O-ring in the half round groove in the bottom plate.

Insert the sample to be desiccated.

Place the clear desiccator top onto the base, making sure that the bottom edge comes into contact with the O-ring. Press down until the flange snaps into the two locks.

To remove the top, turn the valve to shut off the vacuum and then turn to vacuum release position. Press back one of the snap locks and raise the top section.




OTHER FINE SCIENCEWARE® PRODUCTS FOR YOUR LABORATORY

Demand More From Your Desiccator



Secador™ 1.0
Volume .75 cubic feet

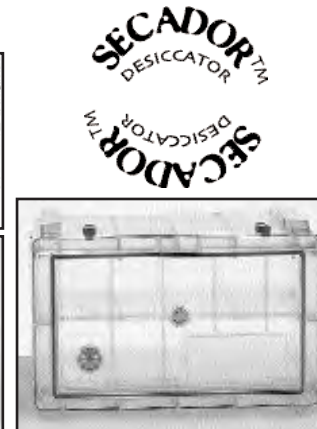

Secador™ 4.0 Horizontal & Vertical
Volume 2.0 cubic feet

Select the Secador™ Desiccator Cabinet that suits your lab!

- Patented sealed construction – Airtight, dust and moisture free storage
- Heavy duty copolyester composition builds rugged cabinet with long service life
- Imbedded UV protection blocks 99% of light from cabinet and contents
- Superior Chemical Resistance - Prevent staining and crazing
- Built in Hygrometer for easy RH monitoring
- Stackable – Saves valuable bench space
- Use with any Scienceware® desiccant

Description	Catalog Number
Secador™ 1.0 Desiccator Cabinet	42071-0000
Secador™ 4.0 Desiccator Horizontal Cabinet	42074-0000
Secador™ 4.0 Desiccator Vertical Cabinet	42074-1000

Bel-Art Products

661 Route 23 South, Wayne, NJ 07470-6814 USA
TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • www.belart.com

94109-9001
Scienceware® is a reg. TM of Bel-Art Products • No-Wire™ ...trademark of Bel-Art Products
Bel-Art Products assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products. All such advice is given and accepted at the buyer's risk.

10/02



Bel-Art Products

661 Route 23 South, Wayne, NJ 07470-6814 USA
TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • www.belart.com