


INSTRUCTIONS



Secador® Auto-Desiccator Cabinet Replacement Desiccant

Catalog No.
42074-0020

1. Disconnect the electronic desiccating unit located on the door of the Secador® Desiccator cabinet by pulling out the electrical connector on the lower edge.
2. Release the electronic desiccating unit from the door by unscrewing the two holding screws on the ends a few turns. Do not remove the screws.
3. Do not touch the four window flaps or metal retainers of the unit. Remove the unit and place it on the lab bench with the soft plastic plug upward.
4. Remove the plastic plug and discard.
5. Pour the used desiccant out of the unit without disturbing the flaps or metal retainers.
6. Remove the cap from bottle containing new desiccant and insert the neck of the bottle into the well from which the soft plastic plug was removed. Pour the entire contents of the bottle into the unit.
7. Insert the new replacement plastic plug into the filler hole, being sure that it is completely inserted.
8. Attach the electronic desiccating unit to the Secador® cabinet door. **Be sure that the washers are directly under the screw heads.** Do not overtighten the screws.
9. Reconnect the electrical connector inside the cabinet to the electronic desiccating unit.

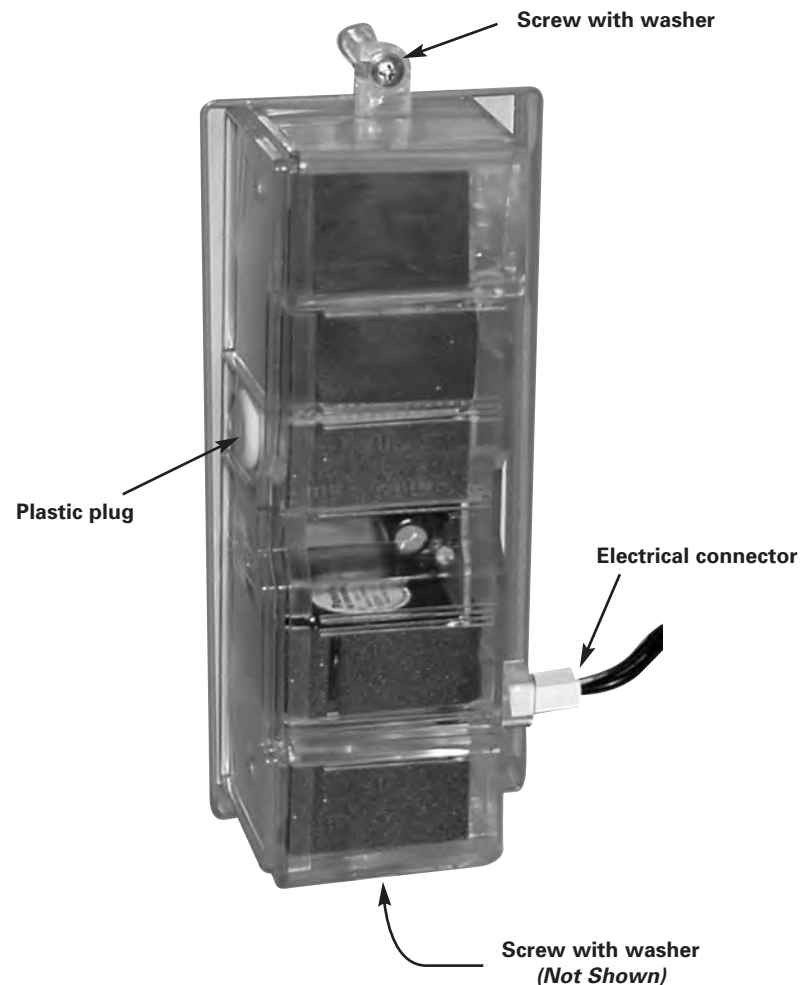

661 Route 23 South, Wayne, NJ 07470-6814 USA
TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • www.belart.com

942070114

05/15

Bel-Art – SP Scienceware assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products. All such advice is given and accepted at the buyer's risk.

© 2015. All Rights Reserved.

Bel-Art® – SP Scienceware® is a division of SP Industries, Inc.