



Floating Bubble Rack
Catalog No. F18875-1600, 1000, 0400



USE AS FOLLOWS:

To Fill:

1. Place the rack on a bench top.
2. Insert the empty tubes.
3. Fill tubes.
4. Close caps securely
5. Lift rack slightly & push down on tubes to lock securely in rack.
6. The rack is now ready for flotation

To Empty:

1. Place filled rack on the bench.
2. Push down the rack to release tubes.
3. Remove tubes.

Let others know what you think. Write a review at www.belart.com

661 Route 23 South, Wayne, NJ 07470

TEL: 973-694-0500 • FAX: 973-694-7199 • www.belart.com

Bel-Art – SP Scienceware assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products. All such advice is given and accepted at the buyer's risk. © 2018. All Rights Reserved.

Bel-Art® – SP Scienceware® is a division of SP Industries, Inc.



999320672 – 09/18



YOU MAY ALSO LIKE

Microcentrifuge Tube Refrigerator Rack
Keep Your Microtubes Cool!



Cat. No. H18901-0000

Cat. No. H18907-0000

- *Two Piece Styrene Tray/Rack*
- *Bottom Section Designed to Hold Water or Ice for Stabilizing Temperatures*
- *Large Capacity Stackable Racks*
- *Available in Two Sizes to Hold a variety of 400µl or 1.5ml tubes.*

For more information, visit www.belart.com