


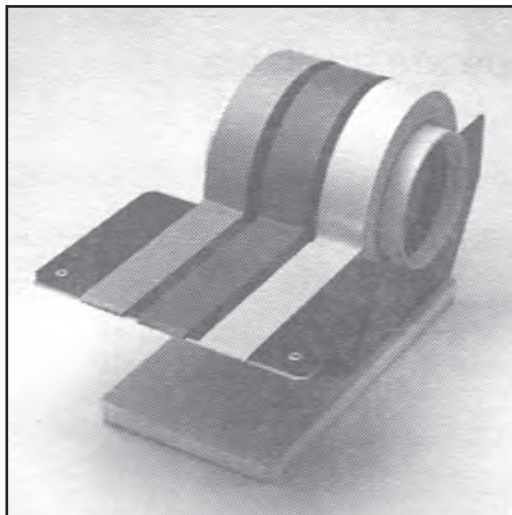

Multiple Roll Tape Dispenser Catalog No. H13468-Series

Remove the white plastic drum to insert rolls of tape. Place spacer rings between rolls of tape to separate them and avoid sticking. Mount all rolls in the same direction so that tape feeds off the top toward the platform with adhesive side down. Install the drum into the frame with slots downward to engage the frame sides. Pull the tape over the platform and write the message, then pull the tape up off the platform and cut on the serrated edge to desired length. To advance the tape, place a finger under the tape between the platform and the drum and lift it until it releases. The dispenser will hold 6 rolls 1/2" wide, 3 rolls 1" wide or 4 rolls 3/4" wide. A combination of widths is also possible.



TO ORDER TAPE: (specify from numbers below)

Catalog No.	W x L / Core	Color
F13483-0050	1/2" x 40yds/3"	White
F13483-0075	3/4" x 40yds/3"	White
F13483-0100	1" x 40yds/3"	White
F13490-0050	1/2" x 60yds/3"	White
F13490-0075	3/4" x 60yds/3"	White
F13490-0100	1" x 60yds/3"	White
F13485-0050	1/2" x 40yds/3"	Yellow
F13485-0075	3/4" x 40yds/3"	Yellow
F13485-0100	1" x 40yds/3"	Yellow
F13484-0050	1/2" x 40yds/3"	Red
F13484-0075	3/4" x 40yds/3"	Red
F13484-0100	1" x 40yds/3"	Red

Catalog No.	W x L / Core	Color
F13486-0050	1/2" x 40yds/3"	Green
F13486-0075	3/4" x 40yds/3"	Green
F13486-0100	1" x 40yds/3"	Green
F13487-0050	1/2" x 40yds/3"	Blue
F13487-0075	3/4" x 40yds/3"	Blue
F13487-0100	1" x 40yds/3"	Blue
F13488-0050	1/2" x 40yds/3"	Orange
F13488-0075	3/4" x 40yds/3"	Orange
F13488-0100	1" x 40yds/3"	Orange


Other Fine Bel-Art - SP Scienceware Products For Your Laboratory

Cover your labels with clear plastic tape

Protective Lab Labeling System covers and protects labels with extra duty clear tape.

Three sizes of tape are available

- F13491-0100 1" x 36 yards
- F13491-0150 1 1/2" x 36 yards
- F13453-0020 2" x 36 yards

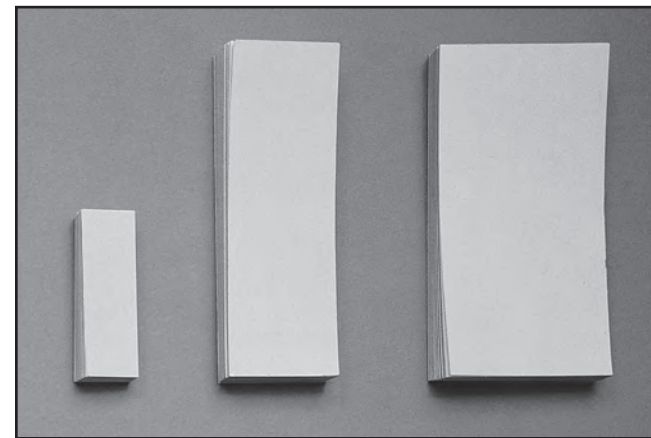

Color coded labels are also available.

White	Yellow	Blue	Green	Dimensions
13455	13456	13457	13458	-
-0005	-0005	-0005	-0005	1/2" x 1-1/2"
-0010	-0010	-0010	-0010	1" x 3"
-0015	-0015	-0015	-0015	1/2" x 3"

661 Route 23 South, Wayne, NJ 07470

TEL: 1-800-423-5278 • FAX: 973-694-7199 • www.belart.com

999320839 - 05/15

Bel-Art - SP Scienceware assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products.

All such advice is given and accepted at the buyer's risk.

© 2015. All Rights Reserved.

Bel-Art® - SP Scienceware® is a division of SP Industries, Inc.

