

I N S T R U C T I O N S



Scienceware® Disposal Carton Covers

Catalog No. F13204-0000 and F13204-0001



Use with corrugated cardboard disposal cartons such as F13205-0001, F24653-0001 and F24657-0001 that have a 30 x 30cm (12 x 12") top opening:

Construct cardboard carton and lid and install plastic liner according to instructions printed on the carton.

Invert the cardboard lid and store on the bottom of the carton.

Place the Disposal Carton Cover over the carton opening. Make sure the lid is oriented so that the side with the release tab is directly over one of the flat edges of the carton.

Gently press the cover down onto the cardboard carton.

Test the spring-actuated pop-up panel by pressing the release tab. Panel should open and remain in open position.

Close the pop-up panel by pushing it down and allowing the release tab to catch.

When disposal carton is full and ready for disposal, remove the cover and replace it with the cardboard lid which was stored at the bottom of the carton. Dispose of carton accordingly.

Disposal Carton Cover may be washed and autoclaved as needed prior to reuse. (see instructions below).

Cleaning Instructions:

IMPORTANT: Before washing, release the panel from its closed position. This releases the spring tension and prevents panel from warping due to exposure to high temperatures.

Can be safely washed with standard laboratory cleaners.

Autoclaving Instructions:

IMPORTANT: Before autoclaving, release the panel from its closed position. This releases the spring tension and prevents panel from warping due to exposure to high temperatures. Recommended autoclave cycle is 121°C at 15 psig (1 bar) for 20 minutes.

Product may be steam autoclaved as needed. It is recommended to wash product before autoclaving to prevent baking contaminants into the surface of the plastic during the autoclave process.

Bel-Art Products

661 Route 23 South, Wayne, NJ 07470 USA

Tel: 1-800-4BELART • Fax: 973-694-7199 • www.belart.com

913192019

08/08

Scienceware® is a registered TM of Bel-Art Products • Bel-Art Products assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products. All such advice is given and accepted at the buyer's risk. ©2008 Bel-Art Products. All Rights Reserved.



Bel-Art Products

661 Route 23 South, Wayne, NJ 07470 USA

Tel: 1-800-4BELART • Fax: 973-694-7199 • www.belart.com