

# I N S T R U C T I O N S



## Benchtop Biohazard Disposal Can Catalog No. H13194-1011

### High-Performance Automation

Traditional bins require you to touch them with your hand or foot to open, which is both inconvenient and unhygienic. This automatic opening bin, based on infrared technology removes this need. Simply hold your hand about 10cm above the infrared panel and the lid will open automatically. Within a few seconds, it will close again automatically.

### Instructions:

- 1. Installation:** Open the battery compartment in the bin and insert batteries (3 AAA - not included).
- 2. Turn on** the power switch located next to the battery inserts. The red indicator light will stay on for 5 seconds, then flash once every 3 seconds to show it is in working order.
- 3. Automatic Opening:** Hold your hand about 10cm above the infrared panel for the lid to open. Move your hand away and lid will close automatically within 5 seconds.
- 4. Manual Opening:** Press the left button on the infrared panel. The lid will open and stay open without closing automatically. Press the right button and the circuit system will return to automatic operation.
- 5. Indicator Lights:** The light in its normal state flashes red, every 3 seconds. When the lid opens, the light shows green. When the lid closes, the light shows red. When the lid is manually opened, it flashes green (once every 3 seconds).
- 6. Batteries:** Batteries will last on average for 6 months, on the basis that the bin is opened/closed 20 times each day. When the batteries need replacement, the indicator light will flash yellow.



Battery Compartment

### Cleaning Instructions:

- To clean, simply wipe with a damp cloth.  
(Do not use running water or heavy liquids, as it will damage electronic components)
- Plastic liner and inside of can may be cleaned using Ethanol (70%) or Bleach  
(Please avoid contact with outside of can or top as finish damage can occur)

### Attention:

- Do not press the lid of the bin when it is opening.
- Remove batteries once flat to avoid seeping liquid damaging the electronic components.
- Avoid direct sunlight to the infrared panel.

Let others know what you think. Write a review at [www.belart.com](http://www.belart.com)



661 Route 23 South, Wayne, NJ 07470

TEL: 973-694-0500 • FAX: 973-694-7199 • [www.belart.com](http://www.belart.com)

913194007 - Rev 2

Bel-Art - SP Scienceware assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products.

All such advice is given and accepted at the buyer's risk.

© 2018. All Rights Reserved.

Bel-Art® - SP Scienceware® is a division of SP Industries, Inc.



### Bag Recommendation:

- Can be used with Benchtop Biohazard Bags
  - F13166-0000 - Pack of 100
  - F13166-0001 - Pack of 1000



## YOU MAY ALSO LIKE

### Biohazard Incinerator Cartons

Safe Disposal, No Need to Transfer Contents



F13205-0001 - Floor model, 6/Pack

### Autoclave Safe Bag Clips - Clavies®

Lets Pressure Out, Keeps Contents In



F13190-0100 - 100/Pack

### Odo-Clave® Autoclave Deodorizer

Neutralize Autoclave Odors  
with a Choice of Four Fragrances



H13198-0000 - Clove, 100/Pack  
H13198-0001 - Cherry, 100/Pack  
H13198-0002 - Pine, 100/Pack  
H13198-0003 - Lemon, 100/Pack  
H13199-0000 - 3x Clove Scent, 100/Pack

### Biohazard Autoclave Gloves - Clavies®

Safely Unload Your Autoclave



H13201-0000 - 33cm (13")  
H13201-0001 - 47cm (18.5")

For more information, visit [www.belart.com](http://www.belart.com)