



## WIRE RACK I.D. TAGS

Catalog No. F13446-0000

### APPLYING & REMOVING THE TAGS:

1. There are two sets of clips on back of each I.D. tag. The larger set are for  $\frac{1}{8}$ " wire and the smaller set are for  $\frac{3}{16}$ " wire. Turn the tag so that the clips you need are adjacent to the top edge of the rack (see photo).
2. Using light pressure, snap the clips onto the top tier of the wire rack.  
**Note:** *only one set of clips can be used to affix the labeling plaque to the wire rack.*
3. To remove, gently pull on tag while holding rack.

### Notes:

- Polypropylene tags can be labeled using dry-erase marker, permanent marker or a pressure sensitive adhesive label.
- Can be autoclaved at 121°C (250°F).

