



I N S T R U C T I O N S



PREPSAFE™ MINI FLOATING MCT RACK Patent Pending

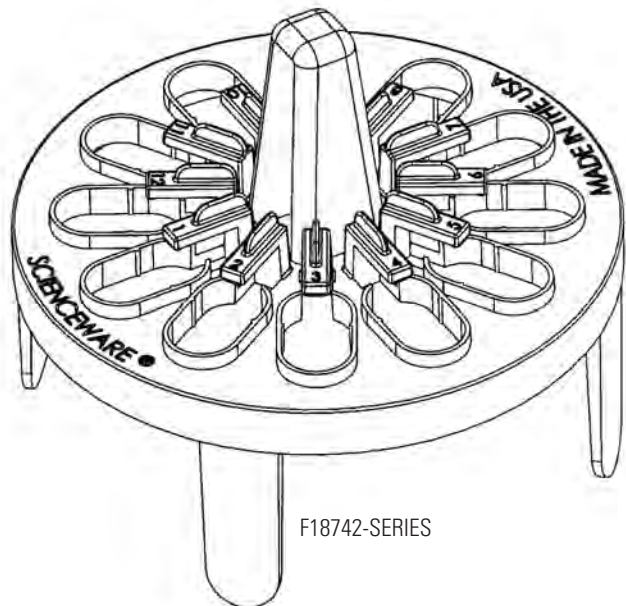

F18742-SERIES

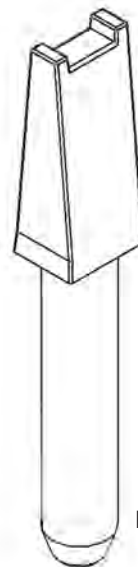

F18742-0140

INSERTING MICROCENTRIFUGE TUBES:

1. Hold rack securely in one hand.
2. Insert a microcentrifuge tube into the desired location.
3. Hold the tube underneath rack and slide it completely under the immobilizing tab until the tube snaps in place securing it in the rack.

REMOVING MICROCENTRIFUGE TUBES:

1. Hold rack securely in one hand.
2. Place your finger underneath the rack behind the tube you wish to remove and slide it out from under the immobilizing tab.

INSERTING VORTEXING ATTACHMENT:

1. Holding the rack in one hand, invert rack so that the legs are pointing upward.
2. Insert the larger end of the vortexing attachment into the hollow space within the handle of the rack until fully engaged.

REMOVING VORTEXING ATTACHMENT:

1. Holding the rack in one hand, invert rack so that the legs are pointing upward.
2. Pull vortexing attachment away from rack.

CATALOG NO.

DESCRIPTION

| | |
|-------------|--|
| F18742-0120 | Clear PrepSafe™ Mini MCT Rack without Vortexing Attachment |
| F18742-0121 | Turquoise PrepSafe™ Mini MCT Rack without Vortexing Attachment |
| F18742-0122 | Fuchsia PrepSafe™ Mini MCT Rack without Vortexing Attachment |
| F18742-0123 | Neon Yellow PrepSafe™ Mini MCT Rack without Vortexing Attachment |
| F18742-0124 | Clear PrepSafe™ Mini MCT Rack with Vortexing Attachment |
| F18742-0125 | Turquoise PrepSafe™ Mini MCT Rack with Vortexing Attachment |
| F18742-0126 | Fuchsia PrepSafe™ Mini MCT Rack with Vortexing Attachment |
| F18742-0127 | Neon Yellow PrepSafe™ Mini MCT Rack with Vortexing Attachment |
| F18742-0140 | Replacement Vortexing Attachment |

