


RETAINS TUBES FOR DECANTING • SUBMERSIBLE

• STACKABLE • AUTOCLAVABLE • SPACE SAVER •

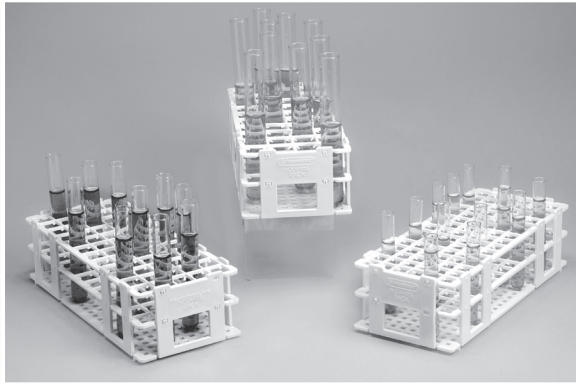

Catalog No.	Holds Tubes	Places	Rows
18749-0000	10-13mm	90	6 x 15
18749-0001	15-16mm	60	5 x 12
18749-0002	18-20mm	40	4 x 10

Assembly instructions on reverse side.



• END PLATES FOR LABELING • EASY TUBE IDENTIFICATION

NO-WIRE™ GRIP RACK



No-Wire™ Grip Test Tube Racks are perfect for every laboratory use, with distinct advantages over other plastic racks. All test tubes are held securely in place, which is especially useful in

decanting the content of all tubes simultaneously. Now, with the addition of this economical new line of No-Wire™ Grip Racks, Bel-Art Products provides you with the widest variety of test tube racks.




661 Route 23 South, Wayne, NJ 07470
TEL: 973-694-0500 • FAX: 973-694-7199 • www.belart.com

I N S T R U C T I O N S



To Assemble:

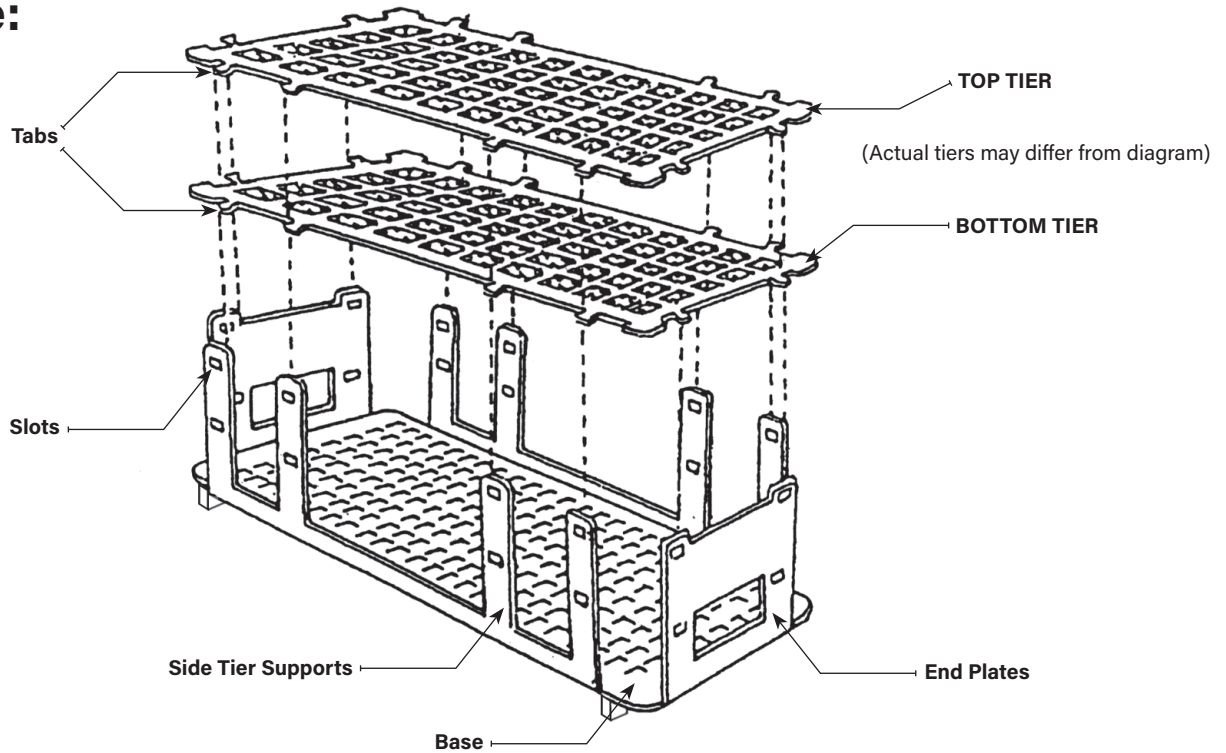

1. Place the base on a flat work surface with the word "BOTTOM" facing down.
2. Bend up one end plate. Snap the bottom tier end tabs into the lower set of slots in the end plate. Be sure the bottom tier imprinted numbers face up.
3. Position the top tier so that imprinted numbers face up and align with the numbers on the bottom tier. Snap the top tier end tabs into the upper set of slots in the end plate.
4. Bend up the opposing end plate and snap the bottom tier and then top tier end tabs into their matching slots in the end plate.

5. Bend up one side tier support and snap the bottom tier and then top tier side tabs into their matching slots in the side tier supports.
6. Bend up the remaining side tier support and snap the bottom tier and then top tier side tabs into their matching slots in the side tier supports.

Submersible and works well in water baths

Autoclavable at 250° F (121°C)

Empty racks can be stacked.

Note: Do not use with liquid nitrogen.

Let others know what you think. Write a product review at www.belart.com

661 Route 23 South, Wayne, NJ 07470

TEL: 973-694-0500 • FAX: 973-694-7199 • www.belart.com

Bel-Art – SP Scienceware assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products. All such advice is given and accepted at the buyer's risk. No-Wire™ is a trademark of Bel-Art – SP Scienceware © 2019. All Rights Reserved.

Bel-Art® – SP Scienceware® is a division of SP Industries, Inc.



918749000 - Rev 2

