




TEST TUBE CAPPER

Catalog No. F18799-0000

Patent Pending

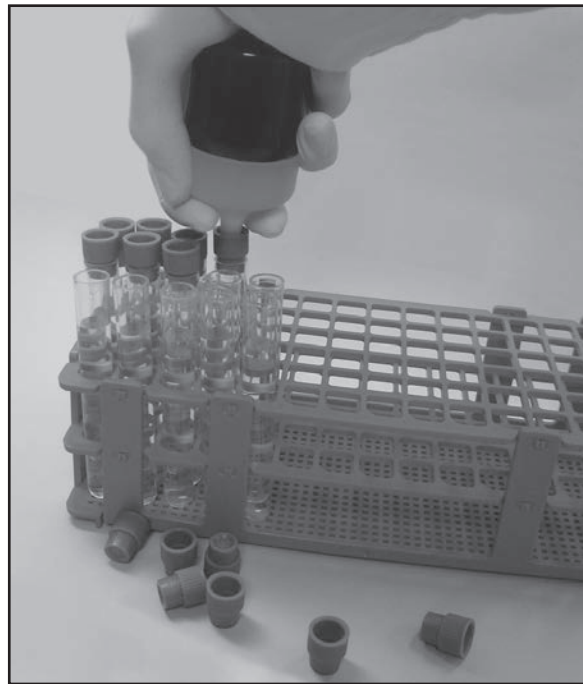

- Eliminate fatigue and soreness of hands and fingers from repetitive task of applying test tube caps
- Fits in the palm of your hand
- Designed for use with recessed center test tube caps

INSTRUCTIONS FOR USE:

1. Grasp capper with the black top resting in the palm of your hand.
2. Place cap on the test tube.
3. Using the Test Tube Capper, gently press cap securely onto the test tube.

661 Route 23 South, Wayne, NJ 07470

TEL: 973-694-0500 • FAX: 973-694-7199 • www.belart.com

Bel-Art – SP Scienceware assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products.

All such advice is given and accepted at the buyer's risk.

© 2016. All Rights Reserved.

Bel-Art® – SP Scienceware® is a division of SP Industries, Inc.



918799002 – 07/16

