

I N S T R U C T I O N S

Maintenance of Scienceware® Eye Wash Bottles

Catalog No. F24850-0000, F24851-0000

1. Wash the Eye Wash Bottle and all accessory parts: cap, tubes, etc. in mild detergent solution.
2. Rinse well to remove all traces of detergent.
3. Under clean conditions, fill Eye Wash Bottle completely.

(We recommend Scienceware® sterile eye wash solution refill # F24879-0032. Please note that once the refill bottle has been opened, the sterility of the solution cannot be assured).

4. Replace top, close cap and return to emergency station.
5. If an eye wash bottle is used, empty it completely and proceed through steps 1-4 before replacing in station.
6. After using an eye wash bottle, the injured individual should see a physician at once for examination and treatment.

To ensure cleanliness, it is advisable to establish a regular eye wash bottle cleaning and eye wash solution replacement procedure. See the refilling guidelines below.



Guidelines for Refilling Scienceware® Eye Wash Bottles

The below recommendations for solution replacement are based on information gathered from sources outside Bel-Art Products to help assist our customers. Bel-Art Products does not guarantee the accuracy of the information gathered.

Rinse solution and frequency of change:

- *Water only – change once a week.*
- *Water and Preservative (varies by manufacturer) - change every two months.*
- *Pre-mixed Eye Wash Solution - change every 4 to 6 months.*
- *Unopened, Sterile Eye Wash Solution (F24879-0032) - until expiration date on bottle.*

Bel-Art Products Sterile Eye Wash solution (F24879-0032) has a shelf life of up to 3 years - refer to expiration date on bottle. However, once the bottle has been opened, the product is no longer sterile. The Sterile Eye Wash Solution is meant for use with Scienceware® Eye Wash Stations (F24878-1032 and F24878-2032).

If using the Sterile Eye Wash Solution (F24879-0032) to refill Scienceware® Emergency Eye Wash Bottles (F24850-0000 and F24851-0000), the solution can no longer be considered sterile once transferred into those bottles. Replace transferred solution every 4 to 6 months.



661 Route 23 South, Wayne, NJ 07470-6814 USA
TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • www.belart.com

Eye Wash Station and Refill



Catalog No. F24868-0000



Catalog No. F24879-0032

Bright, Easy to Locate

Our easily recognized Eye Wash Stations come complete with eye wash bottles, mirror, and instructions for use and service. A bracket holds the bottles securely in place, yet the bottles are easily released when pulled out. Includes station and empty eye wash bottles. We recommend filling with sterile eyewash solution catalog number 24879-0032.

Catalog No.

F24865-0000
F24866-0000
F24868-0000

Bottle Size

Station with one 500ml (16oz) empty eye wash bottle
Station with one 1000ml (32oz) empty eye wash bottle
Station with two 1000ml (32oz) empty eye wash bottles

Recommended Refill

(32oz) sterile eyewash solution refill

F24879-0032



661 Route 23 South, Wayne, NJ 07470-6814 USA
TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • www.belart.com