

I N S T R U C T I O N S



Battery Powered Magnetic Stirrer

F37017-0000



Battery Installation:

1. Separate top half of the stirrer (white) from the bottom half (blue).
2. Install two (2) D size batteries ensuring that both the positive (+) and negative (-) are oriented as shown in the holders.
3. Place top half of the stirrer (white) onto the bottom half (blue), and snap into place by applying slight pressure.

Operation:

1. Ensure that the power switch is in the **OFF** position and that the speed control knob is set at the lowest speed.
2. Place vessel, containing the liquid to be stirred and a Spinbar[®] Magnetic Stirring Bar, on top of the stirring platform.
Note: For optimal performance, do not use Spinbar[®] Magnetic Stirring Bars longer than 1".
3. Center the vessel and the Spinbar[®] Magnetic Stirring Bar on the stirring platform.
4. Turn the power switch to the **ON** position.
5. Adjust stirring speed as needed by using the speed control knob. Rotate the speed control knob clockwise to increase the speed of the stirrer and counterclockwise to reduce the speed.

Caution: When the stirrer is not in use, ensure that the power switch is in the "OFF" position.

Note: For maximum stirring efficiency always specify Spinbar[®] Magnetic Stirring Bars

BEL-ART PRODUCTS

661 Route 23 South, Wayne, NJ 07470-6814 USA
TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • www.belart.com