






Spinbar[®] Capsule Magnetic Stirring Bar

Catalog Number F37152-0000

ASSEMBLY AND USE:

1. Place the Spinbar[®] magnetic stirring bar into the center of the case half.
2. Align the four pins on the top half with four holes in the bottom half and press the two halves of the case together.
3. With the large center hole facing up, place the encased Stirbar into the solution to be stirred.
4. The vessel with Spinbar[®] Capsule Magnetic Stirring Bar can now be placed on a magnetic stirring machine. Be sure to center the stirring bar in the magnetic field of the machine.
5. To clean the unit after use, pull the two halves of the case apart. The case is made of high-impact polystyrene and cannot be autoclaved.

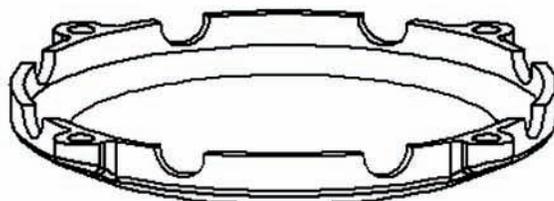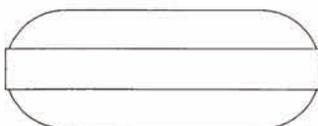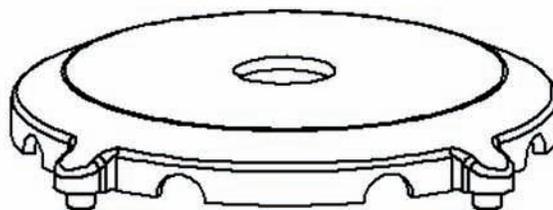

BEL-ART PRODUCTS

Wayne, New Jersey 07470 • Tel: 1-800-4BEL-ART • Fax: 973 694-7199 • www.belart.com