


# INSTRUCTIONS



## Bel-Blotter™

Catalog Number 37876-0002

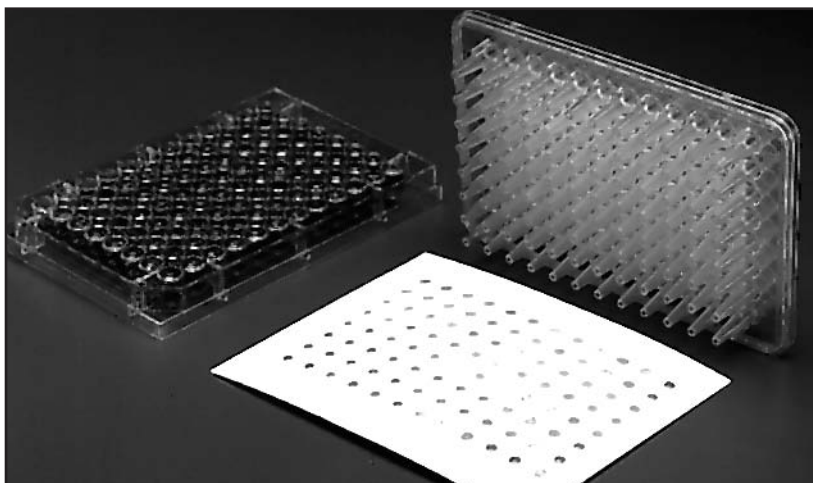

1. The Bel-Blotter, as supplied, is non-sterile, however, it is molded of autoclavable polycarbonate and should be sterilized before use.
2. This replicating tool will do the tedious replicating tasks with little effort and increased productivity.
3. The Bel-Blotter will fit all types of 96 well plates ranging from flat, vee or round bottom plates to the 0.2ml thin walled PCR plates and tubes.
4. By simply placing the Bel-Blotter into the plate, the tips will pick up fluid, up to 10µl per tip.
5. The filled Bel-Blotter is then placed into the receiving medium, whether it be the wells of plates, hybridization membranes or agar.
6. The resulting products of the Bel-Blotter can be used for replication of recombinant DNA libraries, inoculation of filters for colony hybridization, phage typing and several other applications.

## BEL-ART PRODUCTS

661 Route 23 South, Wayne, NJ 07470-6814 USA  
TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • www.belart.com

937876201

02/02

Scienceware® is a reg. TM of Bel-Art Products • Bel-Blotter™ is a Bel-Art Products assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products. All such advice is given and accepted at the buyer's risk.

# OTHER FINE SCIENCEWARE® PRODUCTS FOR YOUR LABORATORY

## Hand Held Electronic Colony Counter

Catalog Number F37862-0000



In a single motion this colony counter marks, counts and confirms with a beep. The lightweight 42.5 gm (1.2oz) pen style colony counter features a four-digit LED display that shows cumulative count. A selector switch allows for counting up or down to correct errors, or to check the count by counting backwards. Powered by a 3V battery. The long life felt-tip marker works on plastic or glass and is oil and water resistant, but removable with ethyl alcohol. Comes with red and black markers. Dimensions: 152.4mm x 19mm x 19mm (6" x 3/4" x 3/4").



## BEL-ART PRODUCTS

661 Route 23 South, Wayne, NJ 07470-6814 USA  
TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • www.belart.com