



Battery Powered Engraving Tool

Engrave all types of materials including glass, plastic, metal and wood



Battery Powered Engraving Tool

- Use to mark and identify test tubes, bottles, racks, equipment, electronics and more
- Engraves most materials including glass, plastic, wood, metal, bone and stone
- Etching is permanent and will not fade with exposure to chemicals, sunlight etc.
- Can be used indoors or outside for field work
- Cordless, uses two 'AA' batteries

Instructions:

The Battery Powered Engraving Tool uses two 'AA' batteries (not included). To insert batteries separate top and bottom sections by gently applying pressure counter clockwise and pulling apart. Place batteries positive side up against spring in lower half and replace top.

Hold the Battery Powered Engraving Tool like you would a pen keeping your thumb over the on/off button. Press the button to start pen and release to stop. To engrave, press head of tool down on the surface at a slight angle and apply light pressure. You may wish to practice on a scrap surface when first using.

To replace tip simply pull out worn tip and push new tip in.

Note: *Always wear protective eyewear when engraving and wash hands after use. Not recommended for engraving electronic media (cd's, dvd's etc.)*

Catalog No. F44150-0005



WAYNE, NJ 07470 USA

TEL: 1-800-423-5278 • FAX: 973-694-7199

www.belart.com