



Cleanware™ Microscope Optics Cleaning Kit Catalog No. F17072-0000

This kit makes it easy to keep optics clear and all microscope parts clean. The kit contains 10 pre-saturated optic and lens cleaning pads, 2 individually wrapped packs of anti-static plastic surface wipes, 2 pre-moistened germicide swabs 6" (152mm) long and complete instructions.

The Value of Regular Maintenance:

Your microscope can give you years of hassle-free use with just a little time spent regularly on cleaning and maintenance. By removing stains from the body and regularly cleaning the optics, you will prevent residue build-up, which causes poor visuals and machine down time.


Cleaning Instructions:

1. Open one lens-cleaning pad and wipe off all optics. No drying necessary.
2. Open resealable pouch and use wiper to clean off entire microscope body.
3. Open one Pre-moistened Swab and clean out all crevices around base, stage and light source.

Cleanware™

Microscope Cleaning Kit

Catalog No. F17072-0000

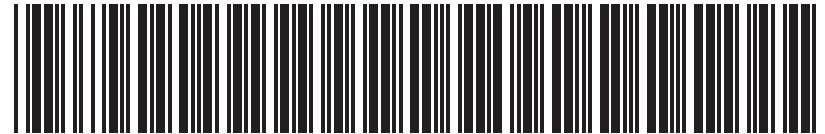

+ H 5 8 8 1 7 0 7 2 0 0 0 0 1

**Everything needed to clean all parts of your microscope
and keep it in optimal working condition**

CONTENTS

- 10 pre-saturated optic and lens cleaning pads.
- 2 pre-moistened germicide swabs.
- 2 anti-static plastic surface wipes.

661 Route 23 South, Wayne, NJ 07470



TEL: 973-694-0500 • FAX: 973-694-7199 • www.belart.com

999320810 - 04/16

Bel-Art - SP Scienceware assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products.

All such advice is given and accepted at the buyer's risk.

© 2016. All Rights Reserved.

Bel-Art® - SP Scienceware® is a division of SP Industries, Inc.



Cleanware™

OPTICAL CLEANER

F17072-0010



+ H 5 8 8 1 7 0 7 2 0 0 1 0 1

Soft, lint free, non-abrasive, pre-moistened cleaning pads for eyeglasses, copier screens, camera lenses, microscopes, telescopes, binoculars, optical scanners viewers or any delicate surface that needs cleaning. Pack includes 10 pads 5" x 8" (127 x 203mm) folded into a 2" x 3" (51 x 76mm) sealed foil pouch.



Wayne, New Jersey 07470 USA